

GEMÜ



**B600 Tank Valves
in Stainless Steel**



Leading the world in pharmaceutical and biotechnology industry sterilisation processes

GEMÜ is one of the leading manufacturers of valves, measurement and control systems for sterile applications in the pharmaceutical and biotechnology industries. This position is based on GEMÜ's comprehensive investments in application-oriented research & development, amounting to more than 5% of the company's turnover. The versatile product range is supplemented with a wide range of advisory services provided by industry specialists and application experts.

Customized solutions for your project business

GEMÜ provides the optimal solution from a single source. As a system supplier of isolation, actuator and control technology, we can respond very flexibly to your individual project-specific needs. Our worldwide sales network provides fast reaction times, customer oriented service and a committed project management team.



Table of contents

Description of use	4	Sterile sampling	
GEMÜ B600 tank valves - Overview of applications	5	Examples with welded on diaphragm valve	20
GEMÜ B600 tank valves - Standard version	6	Sterile take-off/sampling	
Simple draining/tank bottom valve	7	Fixing via clamp connection	21
CIP/SIP of the tank outlet		Tank bottom valve body	
Examples with welded on diaphragm valve	8	with detachable connection	22
CIP/SIP of the tank outlet		Tank bottom	
Examples with integrated diaphragm valve	9	with integrated agitator connection	23
CIP/SIP of the tank		Special applications	24
Examples with welded on diaphragm valve	10	Installation examples	25
CIP/SIP of the tank		Butt weld connections / Surface finish	26 - 28
Examples with integrated diaphragm valve	11	Clamp bodies	29
Tank bottom outlet/venting valve		Materials and certificates	30
with vertical outlet spigot	12	Certificates and approvals	31
Venting valve		The original GEMÜ seal system	32 - 35
Examples with welded on diaphragm valve	13	GEMÜ B600 tank valves - Selection of operators	36 - 38
Tank outlet as ring main		Positioners and process controllers	39
with two outlet spigots	14	Combi switchboxes and electrical position indicators	
Sterile sampling		for pneumatically operated linear valves	40
from a tank	15 - 16	GEMÜ B600 tank valves	
Sterile sampling		Specification	41 - 42
Example with integrated diaphragm valve	17	Worldwide presence	43
Sterile sampling			
from a tank	18 - 19		

Description of use

Today tank valves are available in a large number of versions. They can be installed or welded into the tank cover, tank wall or tank bottom. Their main functions are for filling, sampling and draining the tank contents. Sometimes these functions are combined in one valve for reasons of process safety and sometimes even extra functions are added such as integrated CIP/SIP connections.

All these GEMÜ valves have one thing in common. They are diaphragm valves whose sealing weir is as close as possible to the tank wall to avoid deadlegs in the tank. The internal of the tank bottom valve body has a specially designed cavity which promotes optimum draining of the tank contents and improves their cleanability and sterilisation. Only this cavity ensures optimised draining of the tank.



Features

- Being very compact, the tank valves are ideal for applications where space is at a premium
- Minimal deadlegs and optimized draining capabilities
- The valve body is machined from a single piece of block material (monoblock – no welded components)
- CIP/SIP capable and sterilisable
- The internal surface contour of the valve body is available mechanically and/or electropolished down to Ra 0.25 µm
- Optimised flow geometry
- Pipe connections such as butt weld spigots, clamps and threaded connections are available in accordance with industrial standards
- Valve body materials are 1.4435/316L, other alloys are available to customer specification or test requirements
- Welding into the tank bottom is simplified by a welding neck (standard 6 mm)
- Available with manual, pneumatic or motorized operators from the GEMÜ modular system
- Optical and/or electrical position indicators are available as actuator instrumentation




GEMÜ B600 tank valves

Overview of applications

CIP/SIP of the tank outlet

Tank bottom valve with welded on or integrated diaphragm valve. The welded on diaphragm valve is turned for optimized draining, if technically possible.


For design examples see pages 8-9.

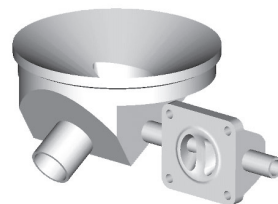

CIP/SIP of the tank or direct sampling from the tank

Tank bottom valve with welded on or integrated diaphragm valve. The welded on diaphragm valve is turned for optimized draining, if technically possible.


For design examples see pages 10-11.

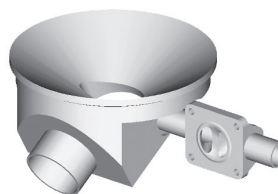

Venting valve

Tank valve with vertical outlet spigot. Available with additional welded on or integrated diaphragm valve depending on the version. The welded on diaphragm valve is turned for optimized draining, if technically possible.


For design examples see pages 12-13.

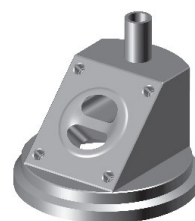

Tank outlet as ring main

Tank bottom valve with two outlet spigots for simple integration into a ring main.


For design examples see page 14.

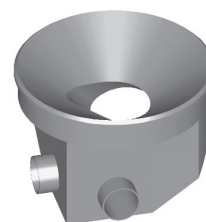

Sterile sampling


Tank valve with integrated sampling option. Also available with integrated CIP/SIP function.

For design examples see pages 15-21.

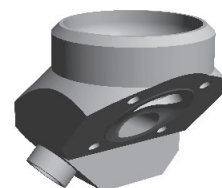

GEMÜ B600 tank valves


Standard version

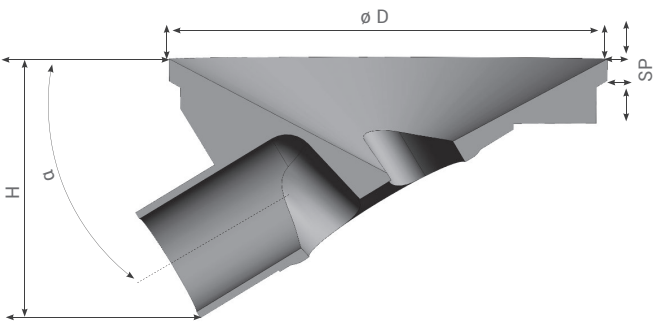

SP = welding neck thickness
 $\varnothing D$ = welding neck diameter
 α = discharge angle
H = valve body height

Diaphragm size	D	SP	H	α
8	$\varnothing 50$	6	29	30°
10	$\varnothing 85$	6	52	25°
25	$\varnothing 120$	6	70	31°
40	$\varnothing 160$	6	95	30°
50	$\varnothing 180$	6	105	33°
80	$\varnothing 250$	6	145	33°
100	$\varnothing 320$	10	198	36°



All dimensions in mm

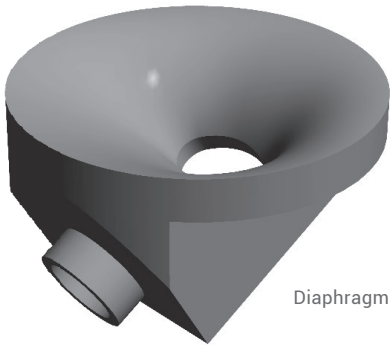


Simple draining/tank bottom valve

Version: Forged design
Application: Pump sump draining

GEMÜ B600 02-01.SK

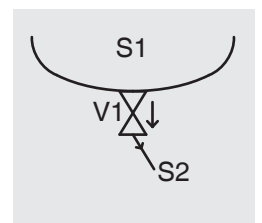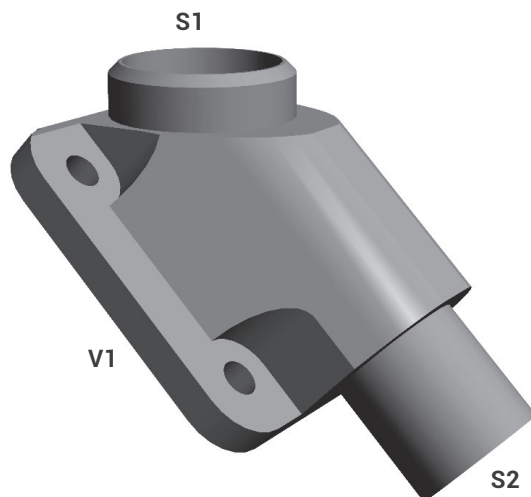

Sectional illustration



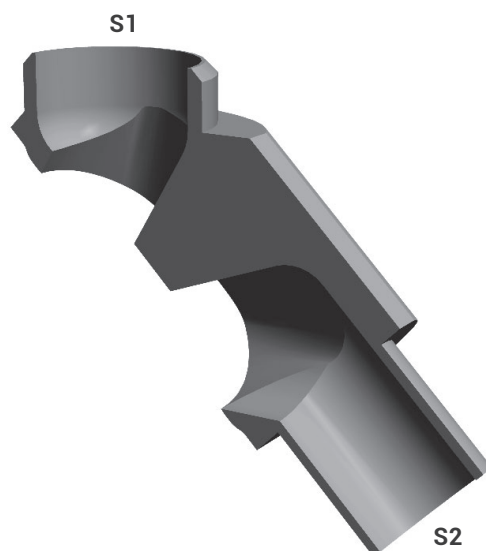


CIP/SIP of the tank outlet




Examples with welded on diaphragm valve

Application: CIP/SIP of the tank outlet
Installation: Tank bottom




GEMÜ B600 03-02.A: Diaphragm valve right, orientation to front

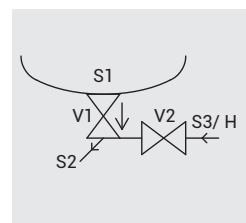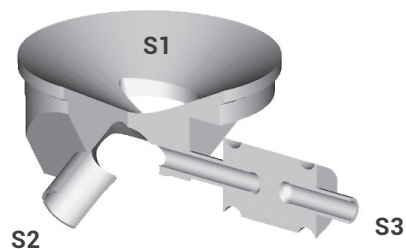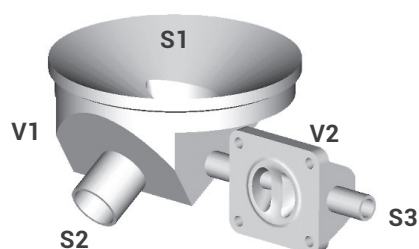




GEMÜ B600 03-02.B: Diaphragm valve left, orientation to front

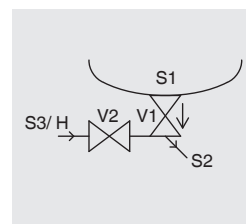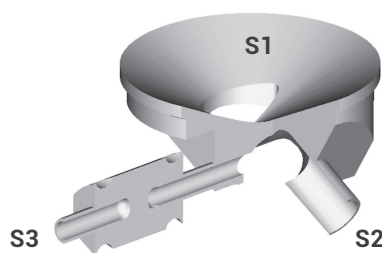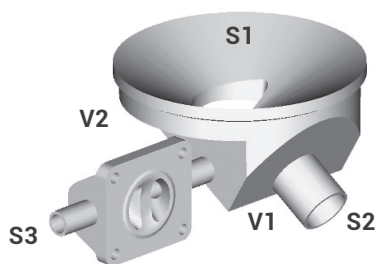

GEMÜ B600 03-02.C: Diaphragm valve right, orientation to back

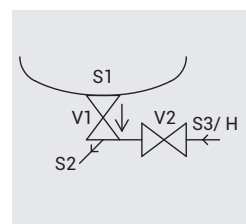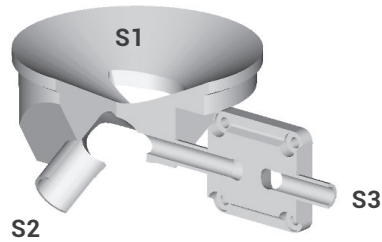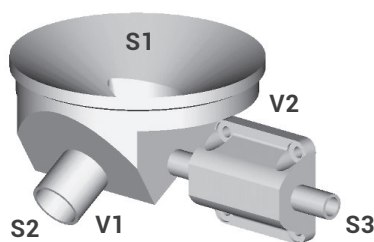

GEMÜ B600 03-02.D: Diaphragm valve left, orientation to back

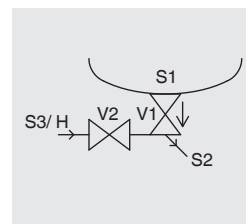
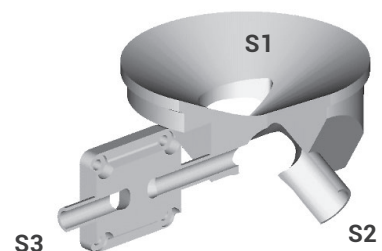
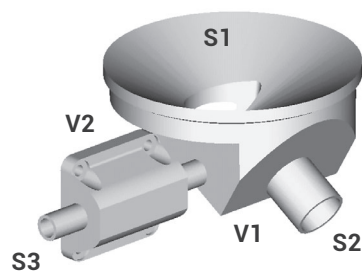


CIP/SIP of the tank outlet




Examples with integrated diaphragm valve

Application: CIP/SIP of the tank outlet
Installation: Tank bottom

GEMÜ B600 03-02.PL

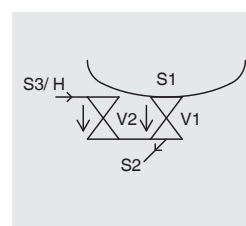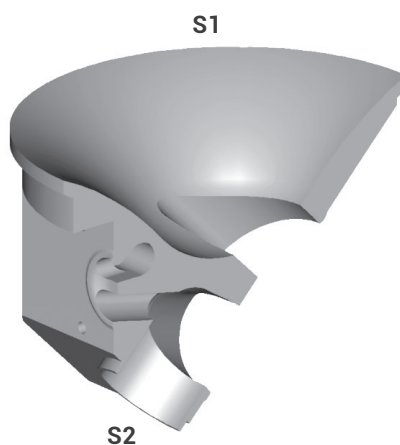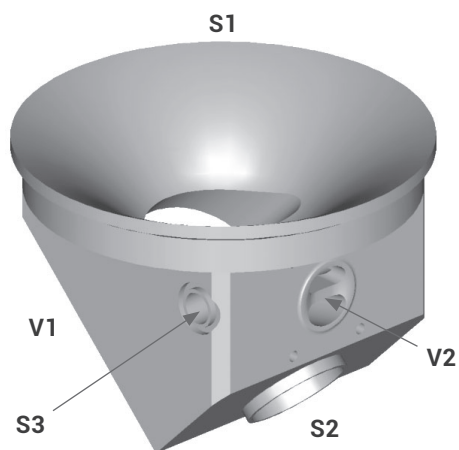


GEMÜ B600 03-02.PR

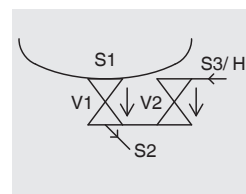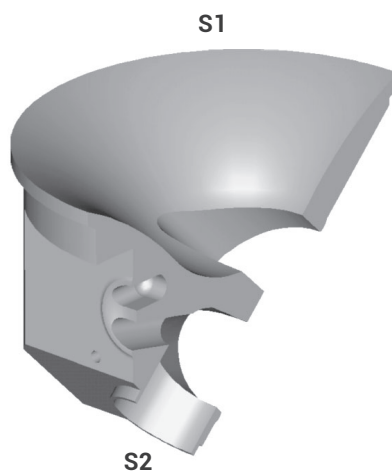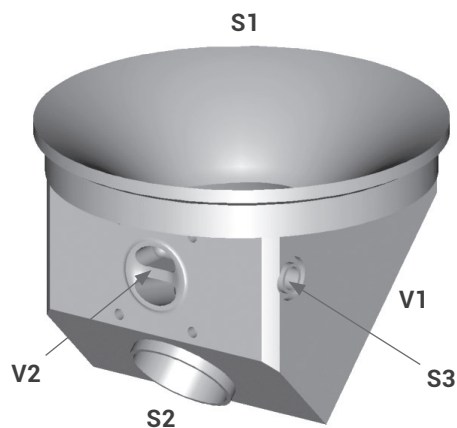

Design for large nominal sizes:

GEMÜ B600 03-02.RR


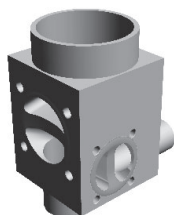


CIP/SIP of the tank



Examples with welded on diaphragm valve

Application: CIP/SIP of the tank or direct sampling from the tank
 Installation: Tank bottom



GEMÜ B600 03-02.K: Diaphragm valve right, orientation to front

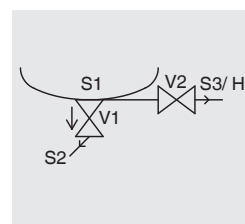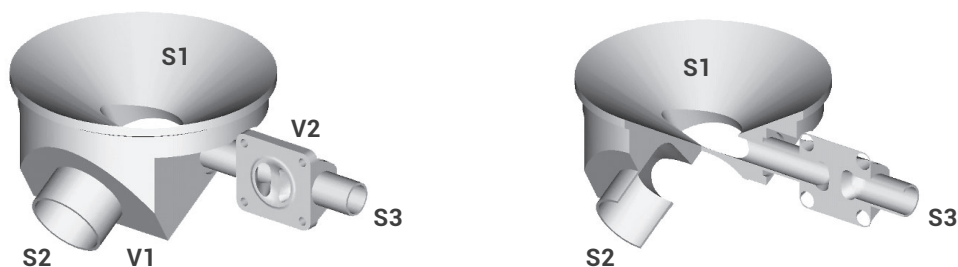



GEMÜ B600 03-02.L: Diaphragm valve left, orientation to front

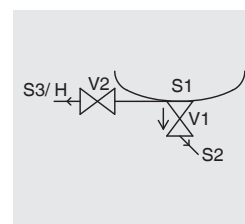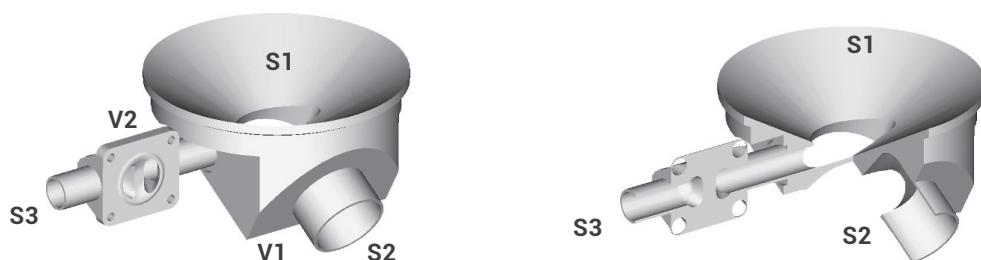

GEMÜ B600 03-02.M: Diaphragm valve right, orientation to back

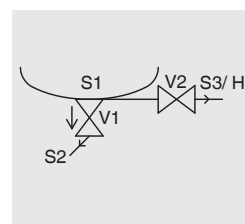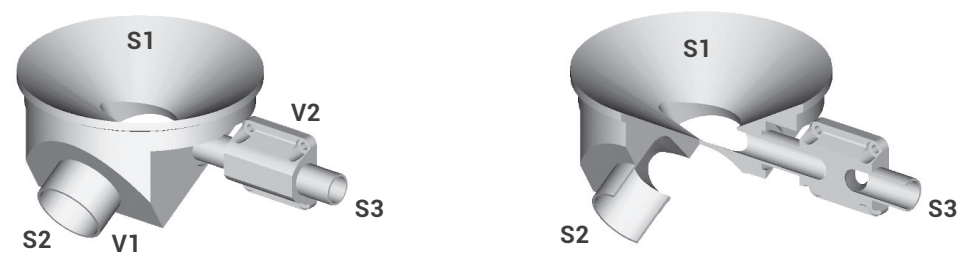

GEMÜ B600 03-02.N: Diaphragm valve left, orientation to back

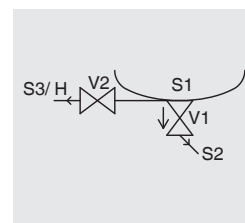
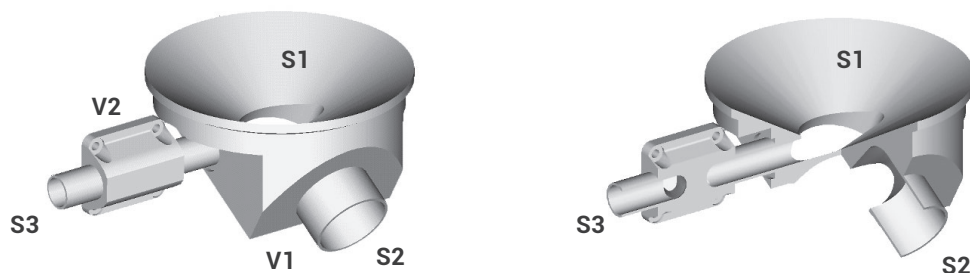


CIP/SIP of the tank




Examples with integrated diaphragm valve

Application: CIP/SIP of the tank or direct sampling from the tank
 Installation: Tank bottom

GEMÜ B600 03-02.SL

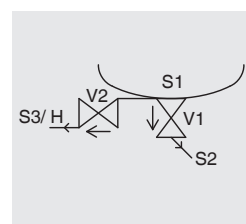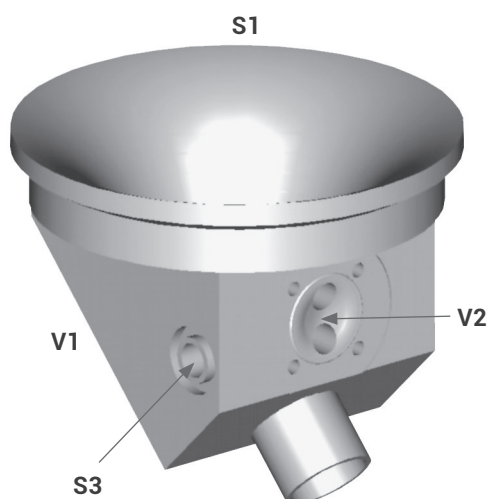



GEMÜ B600 03-02.SR

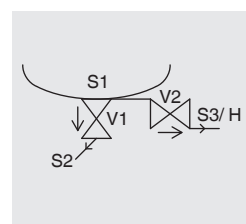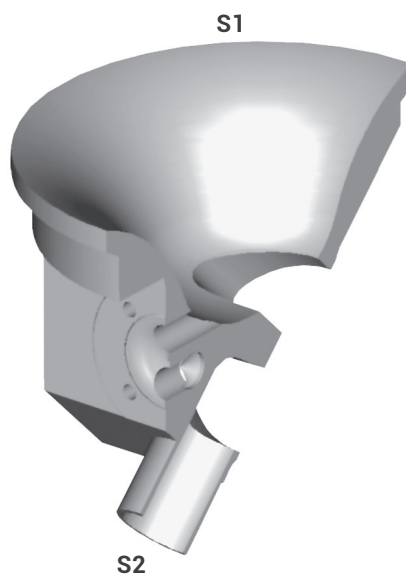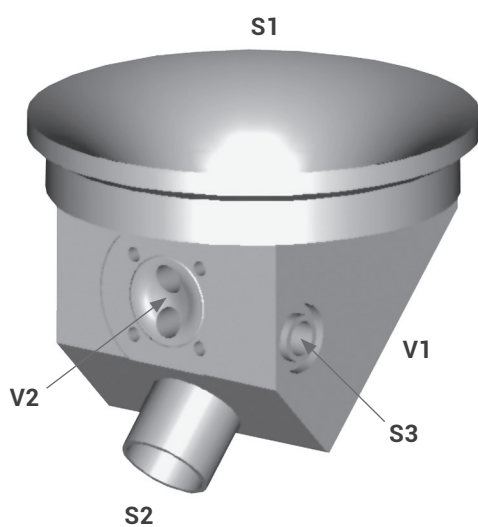


Tank bottom outlet/venting valve with vertical outlet spigot

Application: Outlet valve at the bottom or venting valve in the tank cover
Installation: Tank bottom, tank cover

GEMÜ B600 02-01.H

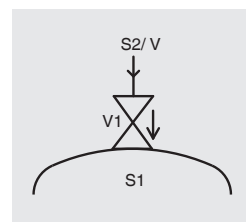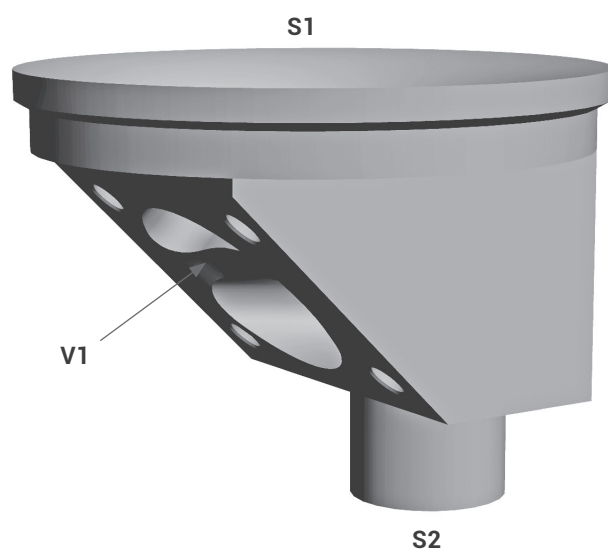

Sectional illustration

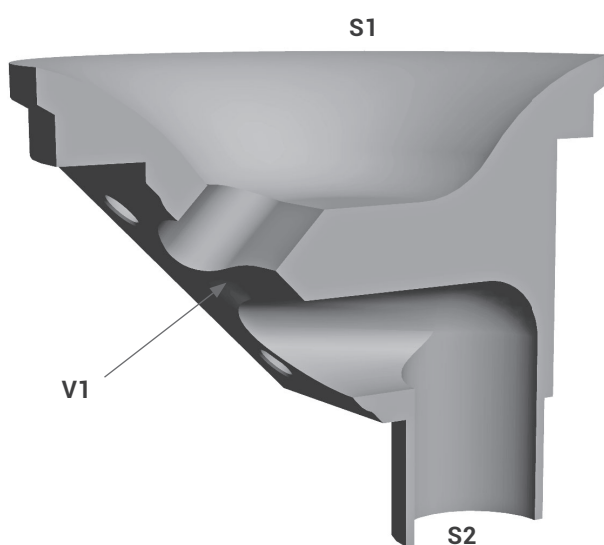

Alternative Einbaulage:





Venting valve

Examples with welded on diaphragm valve


Application: Venting valve in the tank cover with integrated condensate drain or CIP/SIP connection or CIP/SIP of the tank outlet

Installation: Tank cover, tank bottom



GEMÜ B600 03-02.HA: with vertical outlet spigot

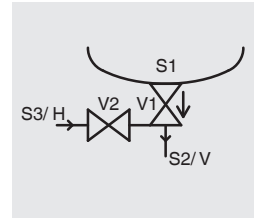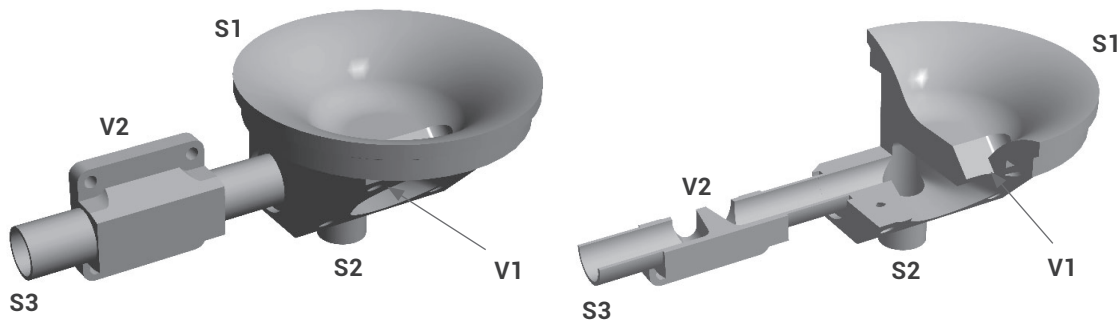


Alternative installation position:

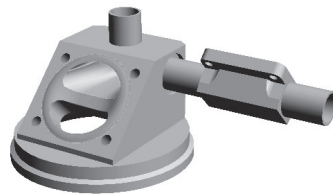

GEMÜ B600 03-02.HB: with vertical outlet spigot

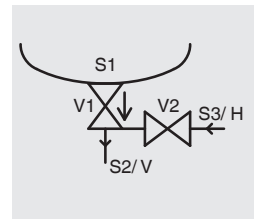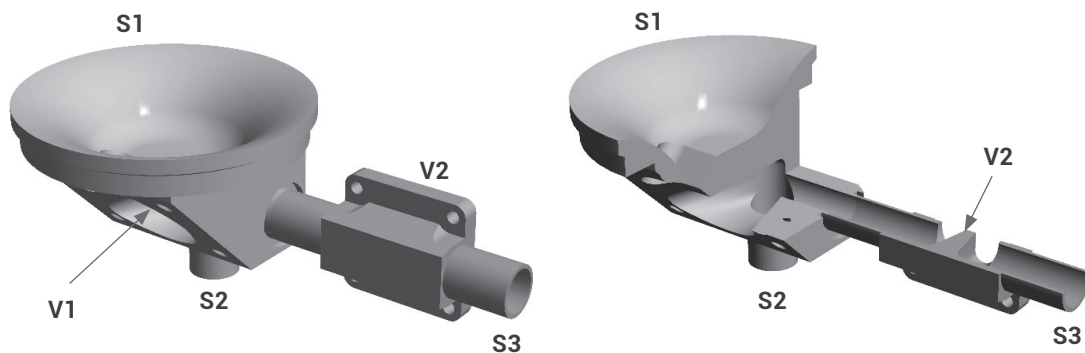




Alternative installation position:

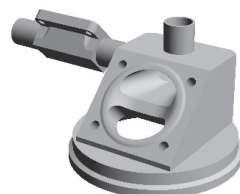




Tank outlet as ring main with two outlet spigots

Application: Tank outlet which can be integrated into ring main
Installation: Tank bottom

GEMÜ B600 03-01.V: Can be integrated into ring main

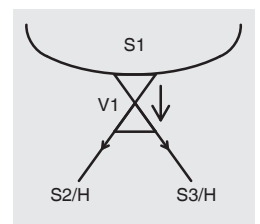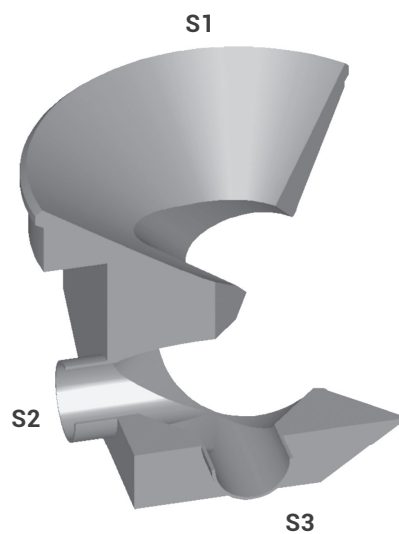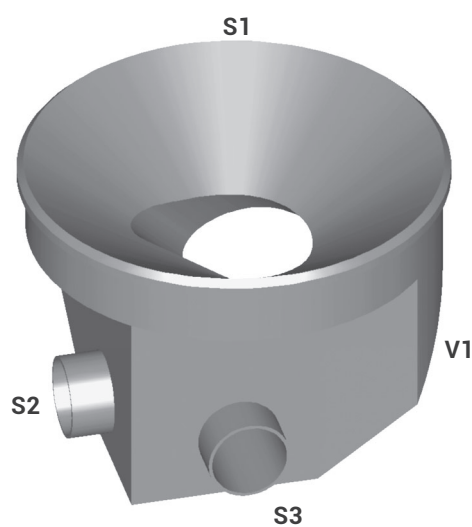




GEMÜ B600 03-01.Y: Ring main directly under the tank

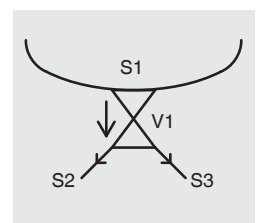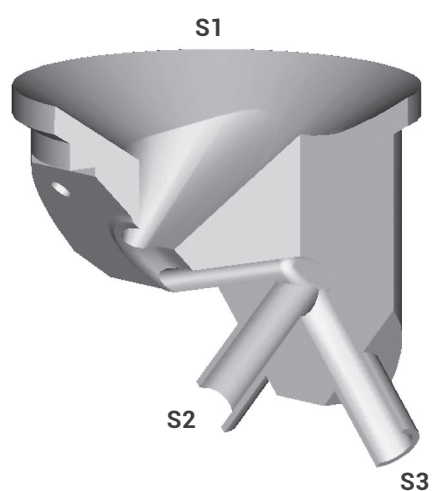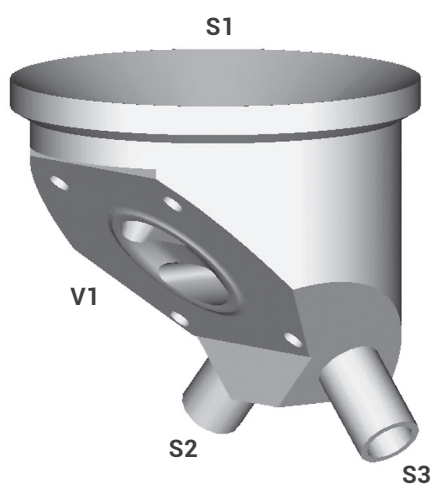




Sterile sampling from a tank

Application: Sampling with integrated flushing connection
Installation: Tank wall
Restriction: Only diaphragm size 8 and GEMÜ 9601, 9602, 9650, 9654 actuators available.
 Max. size of outlets DN 10 acc. to EN ISO 1127.

GEMÜ B600 03-01.Y: Installation suggestion 1

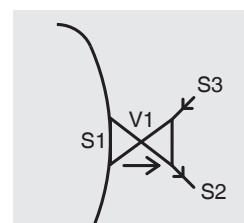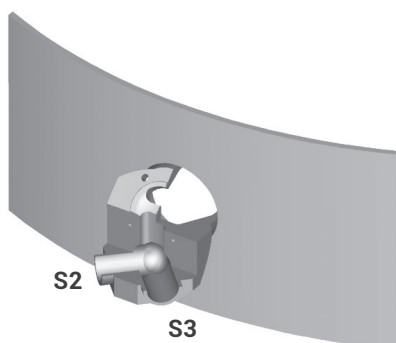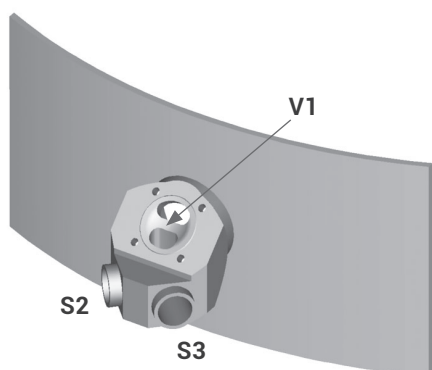




GEMÜ B600 03-01.Y: Installation suggestion 2

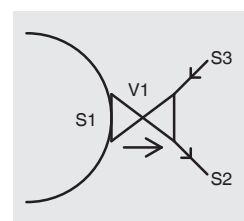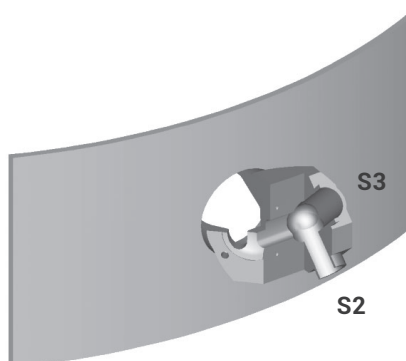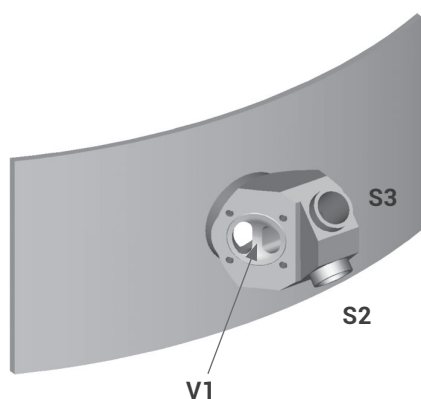




Sterile sampling from a tank

Application: Sampling with integrated flushing connection
Installation: Tank bottom
Restriction: Only diaphragm size 8 and GEMÜ 9601, 9602, 9650, 9654 actuators available.
Max. size of outlets DN 10 acc. to EN ISO 1127.

GEMÜ B600 03-01.Y1: Flushing connection on the right

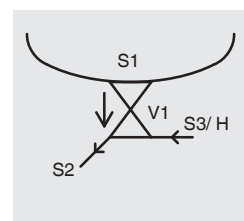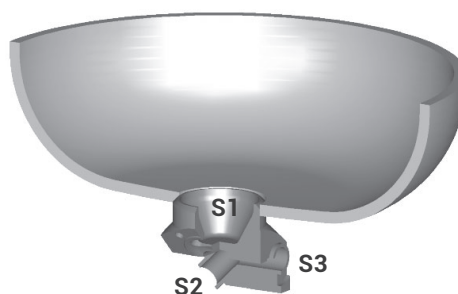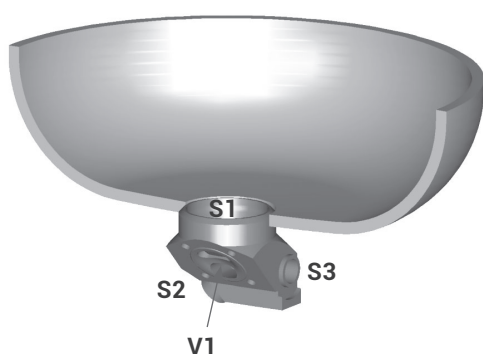

GEMÜ B600 03-01.Y2: Flushing connection on the left

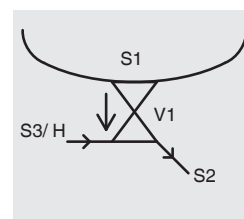
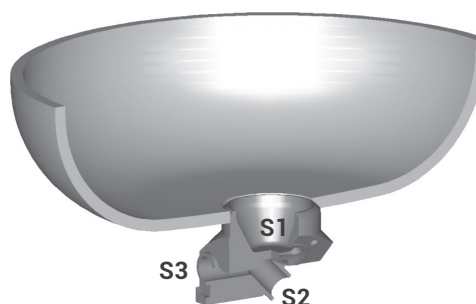
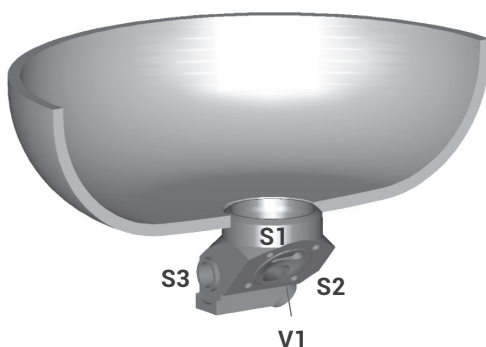


Sterile sampling




Example with integrated diaphragm valve

Application: Sampling with integrated CIP/SIP diaphragm valve
 Installation: Tank wall, tank bottom

GEMÜ B600 03-02.PO

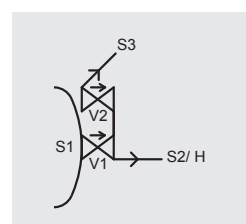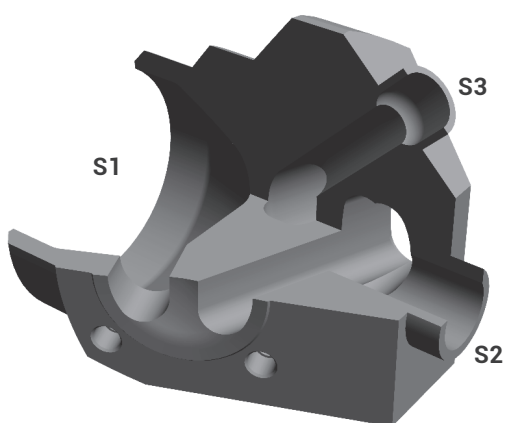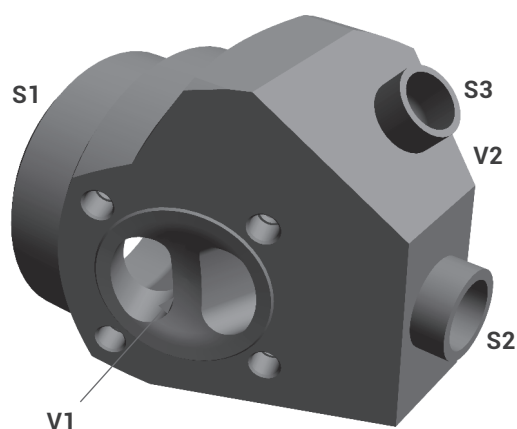


GEMÜ B600 03-02.PU

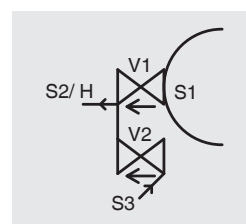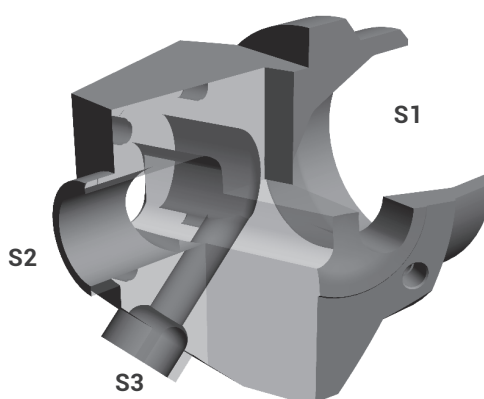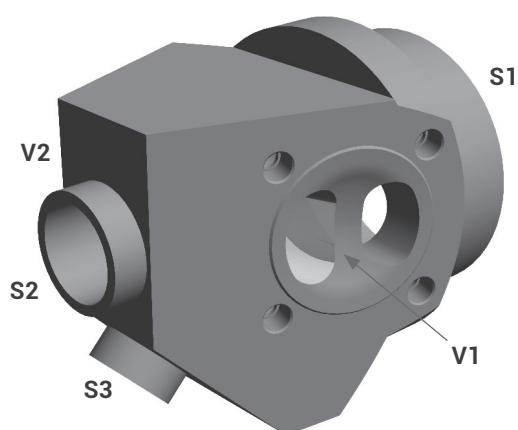


Sterile sampling from a tank

Application: Sampling from a tank
Installation: Tank bottom, tank wall
Restriction: Only diaphragm size 8 and GEMÜ 9601, 9602, 9650, 9654 actuators available.
 Max. size of outlets DN 10 acc. to EN ISO 1127.

GEMÜ B600 02-01.Y1

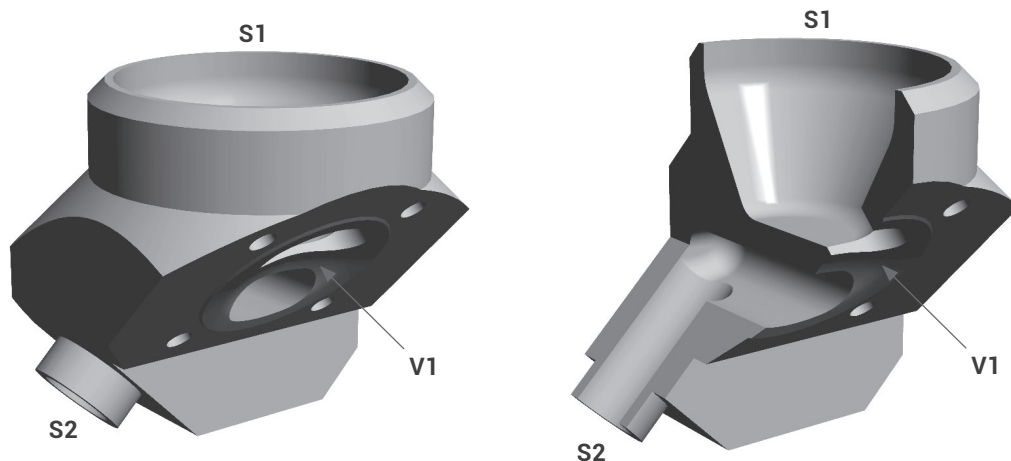



GEMÜ B600 02-01.Y2

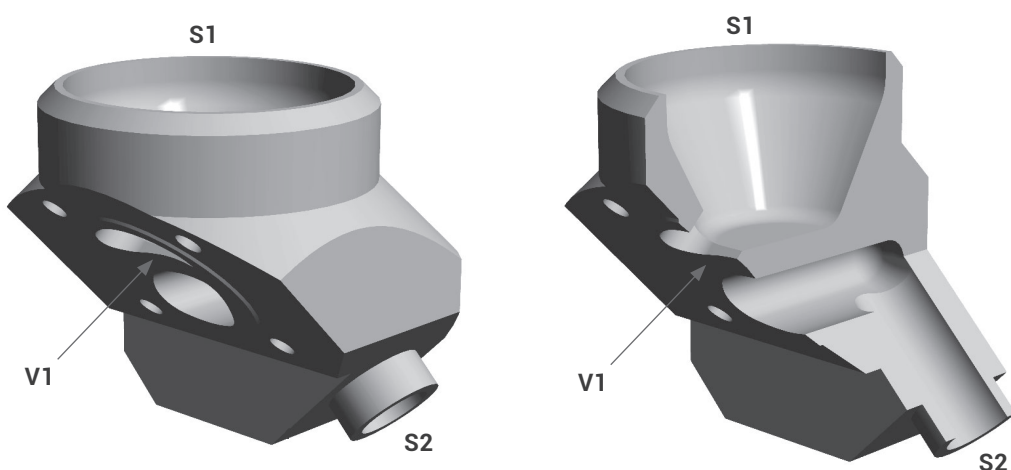


Sterile sampling from a tank

Application: Sampling from a tank
Installation: Tank wall

GEMÜ B600 02-01.F

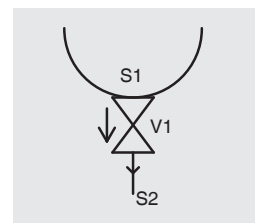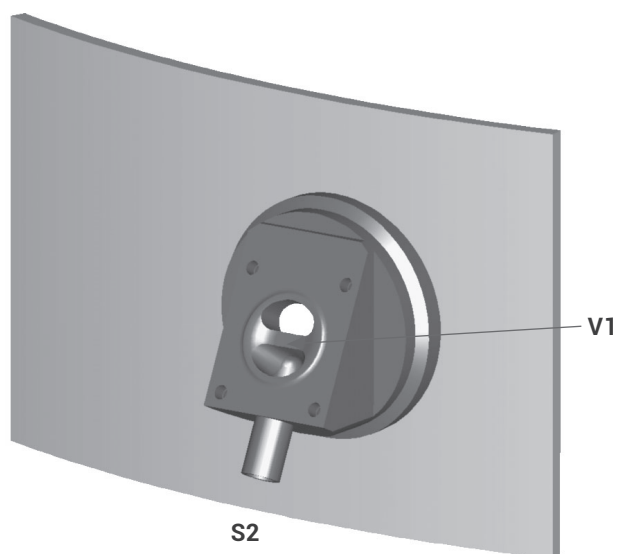

Sectional illustration

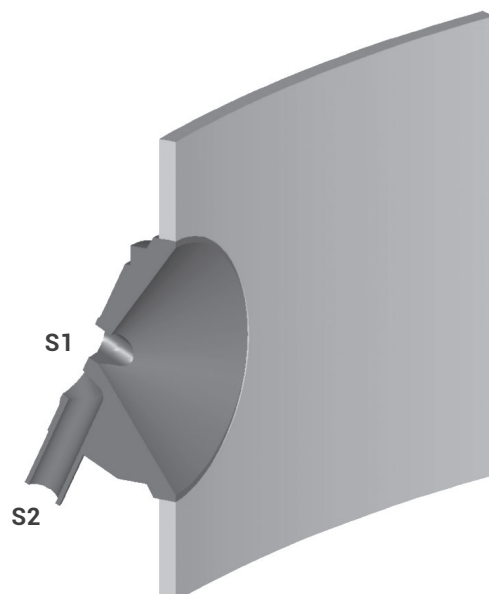


Sterile sampling


Examples with welded on diaphragm valve

Application: Sampling with welded on CIP/SIP diaphragm valve
 Installation: Tank wall

GEMÜ B600 03-02.FL

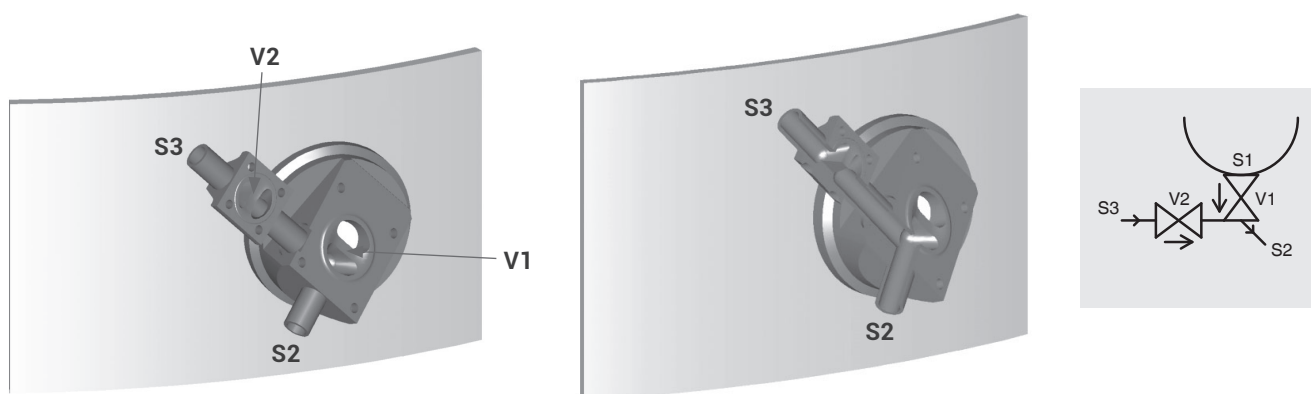

GEMÜ B600 03-02.FR

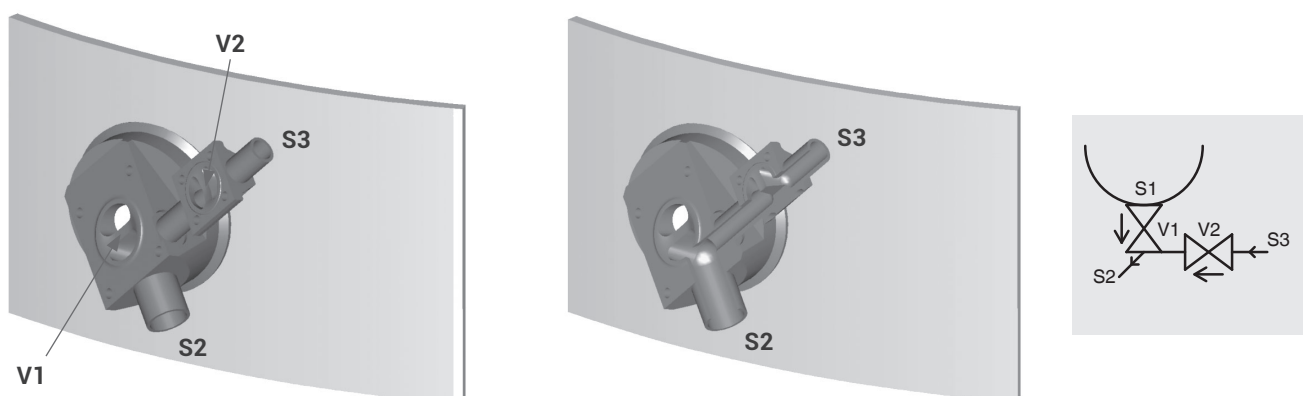


Sterile take-off/sampling


Fixing via clamp connection

Application: Sterile take-off/sampling with integrated clamp connection
Installation: Tank bottom, tank wall
Restriction: Only diaphragm size 8 and GEMÜ 9601, 9602, 9650, 9654 actuators available.
 Max. size of outlets DN 10 acc. to EN ISO 1127.


GEMÜ B600 02-01.Y3

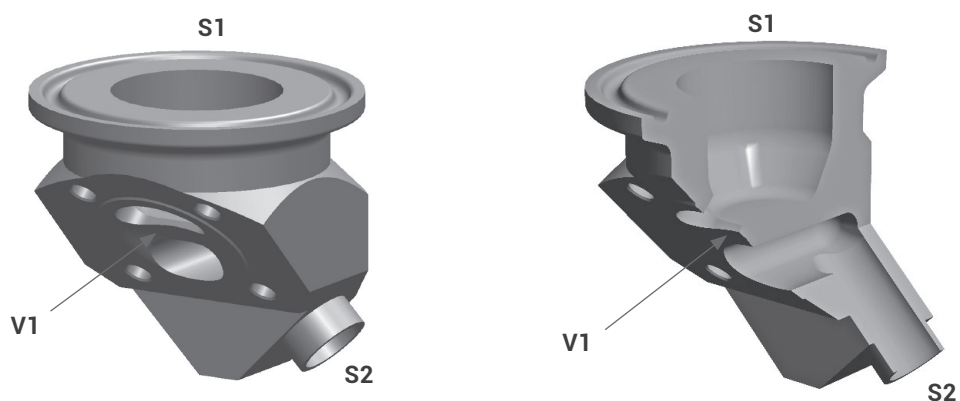


Alternative installation positions:

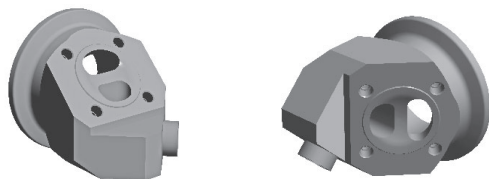

GEMÜ B600 03-01.Y3

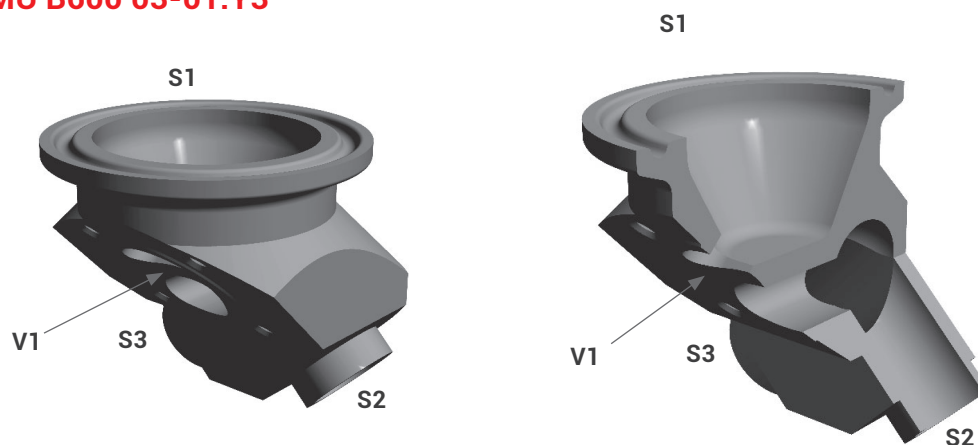



Alternative installation positions:

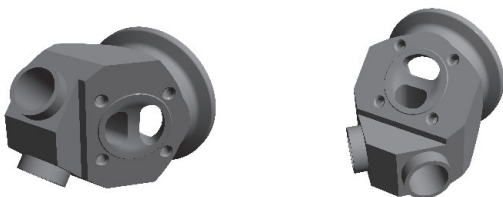


Tank bottom valve body with detachable connection

Application: Sterile take-off/sampling with integrated flanged connection
Installation: Tank bottom

GEMÜ B600 02-01.B

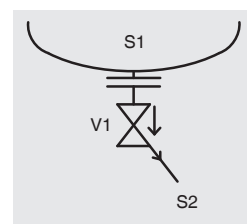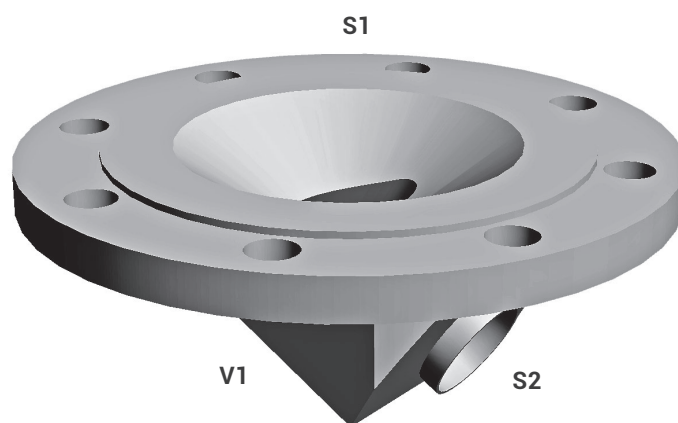

Sectional illustration

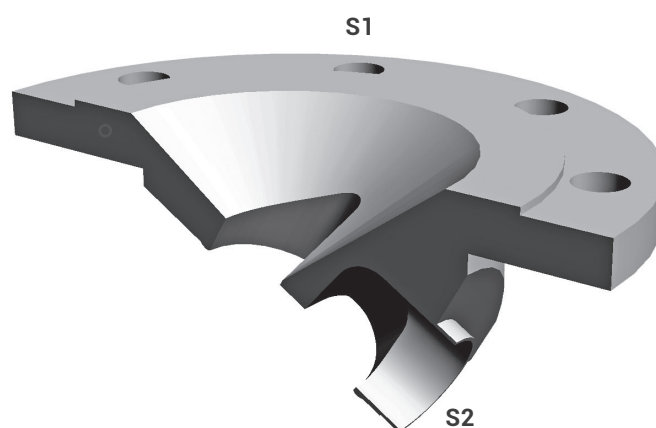




Tank bottom

with integrated agitator connection




Application: Tank bottom outlet with integrated agitator connection (B600 02-01.TB)
or tank bottom outlet with integrated agitator and flushing connection (B600 03-02.TB)

Installation: Tank bottom

GEMÜ B600 02-01.TB

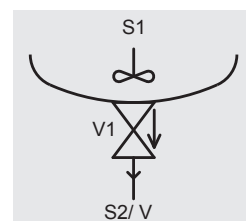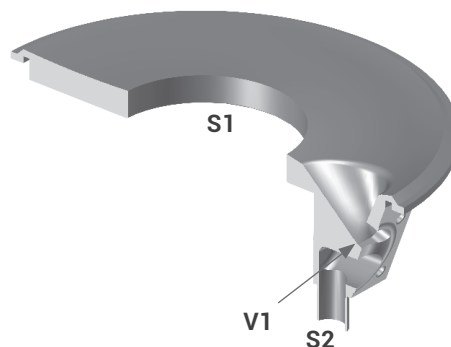
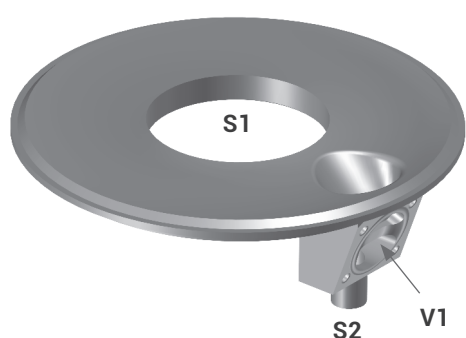


GEMÜ B600 03-02.TB

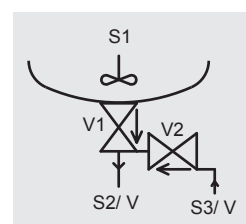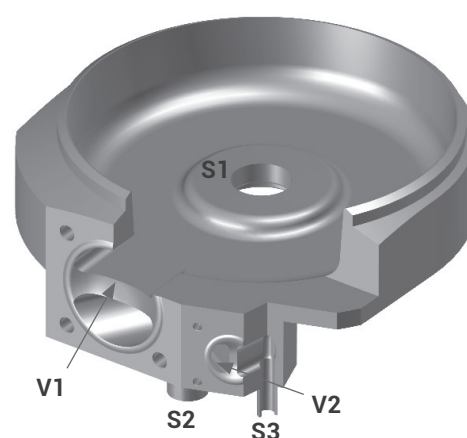
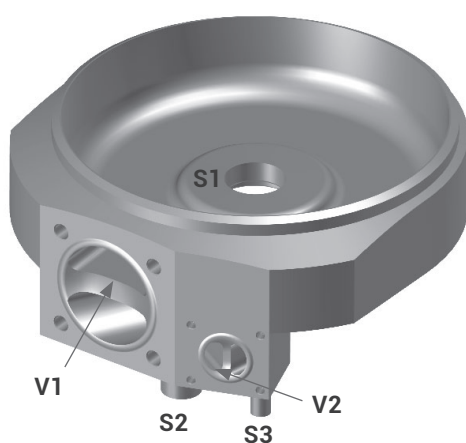


Special applications



GEMÜ B600 07-03.Z

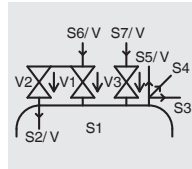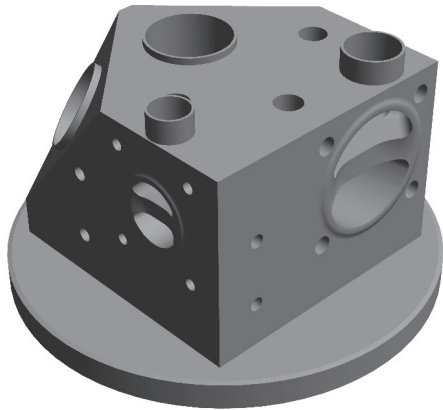



GEMÜ B600 05-06.Z

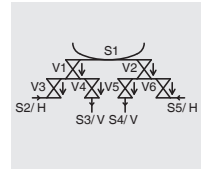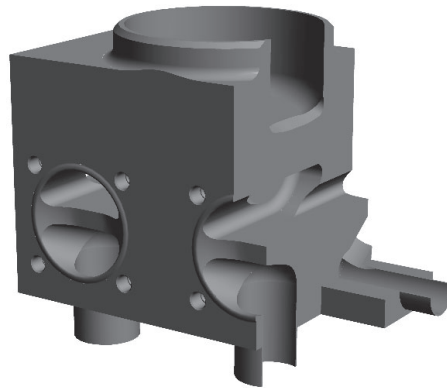



GEMÜ B600 04-04.Z

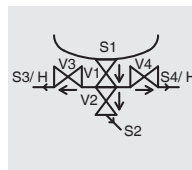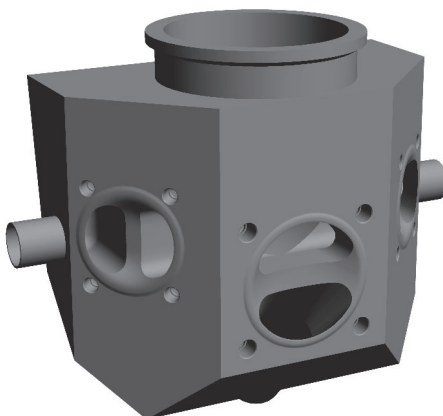

GEMÜ B600 03-03.Z

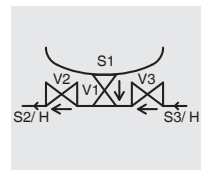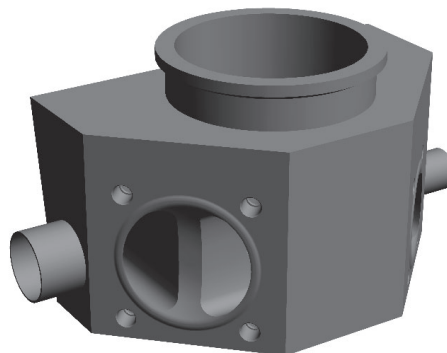

GEMÜ B600 06-04.Z

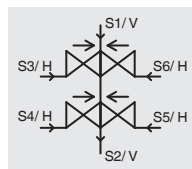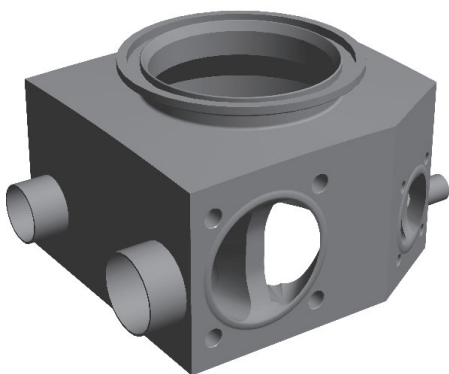

Installation examples



Butt weld connections / Surface finish

Modern, ergonomically shaped workstations and trained polishing staff give us the ability to provide high quality surface finishes. Depending on the required application, surface finishes from Ra 0.8 µm down to 0.25 µm can be achieved by polishing, electro polishing or a special process, we call "elysieren".


Mechanical hand polishing is carried out at our works to ensure our high quality standard.

Valve body surface finish. internal contour			
	Forged body - Codes 40. 42 Block material - Codes 41. 43	Investment casting Codes 32. 34	Code
Ra ≤ 0.8 µm. mechanically polished internal. blasted external	X	X	1502
Ra ≤ 0.8 µm. electropolished internal/external	X	-	1503
Ra ≤ 0.6 µm. mechanically polished internal. blasted external	X ¹	X ¹	1507
Ra ≤ 0.6 µm. electropolished internal/external	X ¹	-	1508
Ra ≤ 0.4 µm. mechanically polished internal. blasted external	X ¹	-	1536
Ra ≤ 0.4 µm. electropolished internal/external	X ¹	-	1537
Ra ≤ 0.25 µm. mechanically polished internal. blasted external	X ¹	-	1527
Ra ≤ 0.25 µm. electropolished internal/external	X ¹	-	1516


Ra acc. to DIN 4768; at defined reference points. Surface finish data refers to media wetted surfaces.

¹ Not possible for connections DN 8 code 55 and 59. DN 4 code 0 and all other connections Ø < 6 mm.

The difference between tube specifications (Example DN 15)

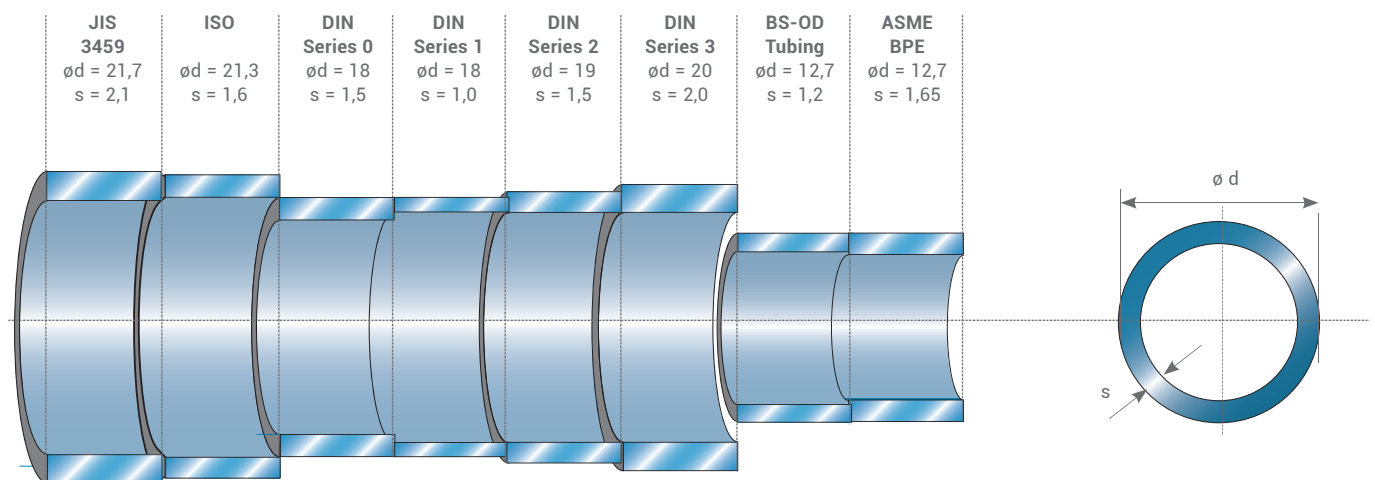

Butt weld connections

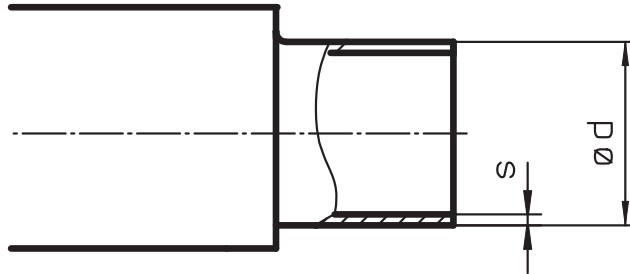

Dimensions in mm			DIN	DIN 11850						DIN 11866						EN ISO 1127	
			Series 0 Code 0	Series 1 Code 16		Series 2 Code 17		Series 3 Code 18		Series A Code 1A		Series B Code 1B		Code 60			
MG	DN	NPS	ød	s	ød	s	ød	s	ød	s	ød	s	ød	s	ød	s	
8	4	-	6	1.0	-	-	-	-	-	-	-	-	-	-	-	-	
	6	-	8	1.0	-	-	-	-	-	-	8	1.0	10.2	1.6	10.2	1.6	
	8	¼"	10	1.0	-	-	-	-	-	-	10	1.0	13.5	1.6	13.5	1.6	
	10	⅜"	-	-	12	1.0	13	1.5	14	2.0	13	1.5	-	-	-	-	
	15	½"	-	-	-	-	-	-	-	-	-	-	-	-	-	-	
10	10	⅜"	-	-	12	1.0	13	1.5	14	2.0	13	1.5	17.2	1.6	17.2	1.6	
	15	½"	18	1.5	18	1.0	19	1.5	20	2.0	19	1.5	21.3	1.6	21.3	1.6	
	20	¾"	-	-	-	-	-	-	-	-	-	-	-	-	-	-	
25	15	½"	18	1.5	18	1.0	19	1.5	20	2.0	19	1.5	21.3	1.6	21.3	1.6	
	20	¾"	22	1.5	22	1.0	23	1.5	24	2.0	23	1.5	26.9	1.6	26.9	1.6	
	25	1"	28	1.5	28	1.0	29	1.5	30	2.0	29	1.5	33.7	2.0	33.7	2.0	
40	32	1 ¼"	34	1.5	34	1.0	35	1.5	36	2.0	35	1.5	42.4	2.0	42.4	2.0	
	40	1 ½"	40	1.5	40	1.0	41	1.5	42	2.0	41	1.5	48.3	2.0	48.3	2.0	
50	50	2"	52	1.5	52	1.0	53	1.5	54	2.0	53	1.5	60.3	2.0	60.3	2.0	
80	65	2 ½"	-	-	-	-	70	2.0	-	-	70	2.0	76.1	2.0	76.1	2.0	
	80	3"	-	-	-	-	85	2.0	-	-	85	2.0	88.9	2.3	88.9	2.3	
100	100	4"	-	-	-	-	104	2.0	-	-	104	2.0	114.3	2.3	114.3	2.3	

MG = diaphragm size

Continued on the next page

Butt weld connections

JIS-G 3447 Code 35					JIS-G 3459 Code 36		SMS 3008 Code 37		BS 4825 Code 55		ASME BPE Code 59		ANSI/ASME B36.19M 10s Code 63		ANSI/ASME B36.19M 40s Code 65	
MG	DN	NPS	ød	s	ød	s	ød	s	ød	s	ød	s	ød	s	ød	s
8	4	-	-	-	-	-	-	-	-	-	-	-	-	-	-	-
	6	-	-	-	10.5	1.20	-	-	-	-	-	-	10.3	1.24	10.3	1.73
	8	1/4"	-	-	13.8	1.65	-	-	6.35	1.2	6.35	0.89	13.7	1.65	13.7	2.24
	10	3/8"	-	-	-	-	-	-	9.53	1.2	9.53	0.89	-	-	-	-
	15	1/2"	-	-	-	-	-	-	12.70	1.2	12.70	1.65	-	-	-	-
10	10	3/8"	-	-	17.3	1.65	-	-	9.53	1.2	9.53	0.89	17.1	1.65	17.1	2.31
	15	1/2"	-	-	21.7	2.10	-	-	12.70	1.2	12.70	1.65	21.3	2.11	21.3	2.77
	20	3/4"	-	-	-	-	-	-	19.05	1.2	19.05	1.65	-	-	-	-
25	15	1/2"	-	-	21.7	2.10	-	-	-	-	-	-	21.3	2.11	21.3	2.77
	20	3/4"	-	-	27.2	2.10	-	-	19.05	1.2	19.05	1.65	26.7	2.11	26.7	2.87
	25	1"	25.4	1.2	34.0	2.80	25.0	1.2	-	-	25.40	1.65	33.4	2.77	33.4	3.38
40	32	1 1/4"	31.8	1.2	42.7	2.80	33.7	1.2	-	-	-	-	42.2	2.77	42.2	3.56
	40	1 1/2"	38.1	1.2	48.6	2.80	38.0	1.2	-	-	38.10	1.65	48.3	2.77	48.3	3.68
50	50	2"	50.8	1.5	60.5	2.80	51.0	1.2	-	-	50.80	1.65	60.3	2.77	60.3	3.91
80	65	2 1/2"	63.5	2.0	76.3	3.00	63.5	1.6	-	-	63.50	1.65	73.0	3.05	73.0	5.16
	80	3"	76.3	2.0	89.1	3.00	76.1	1.6	-	-	76.20	1.65	88.9	3.05	88.9	5.49
100	100	4"	101.6	2.0	114.3	3.00	101.6	2.0	-	-	101.60	2.11	114.3	3.05	114.3	6.02

MG = diaphragm size




Clamp bodies

All clamp connections are machined according to the spigot dimensions e.g. to DIN 11850, EN ISO 1127, SMS 3008 or ASME BPE. We ask our customers to state which version or standard the connections shall comply with.

At GEMÜ all welds are carried out by specially qualified and certified welders utilising state-of-the-art welding technology.

In principle, special connections requested by customers can be provided on GEMÜ M-blocks and it is also possible to have different connections on one body. Valve bodies with fully machined clamp connections are also possible as an option.

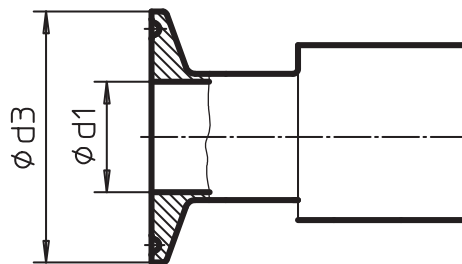

Pipe			Code 59 ASME-BPE				Code 60 EN ISO 1127				Code 59 ASME-BPE				Code 16,17,18 DIN 11850				Code 37 SMS 3008				Code 35 JIS-G3447				Code 36 JIS-G3459			
Clamp connection			Code 80				Code 82				Code 88*				Code 8A				Code 8E				Code 8F				Code 8H			
DN	NPS	MG	H1	ød1	ød3	L	ød1	ød3	L	ød1	ød3	L	ød1	ød3	L	ød1	ød3	L	ød1	ød3	L	ød1	ød3	L	ød1	ød3	L	ød1	ød3	L
8	¼"	8	8	4,57	25	63,5	10,30	25,0	63,5	-	-	-	-	-	-	-	-	-	-	-	-	-	-	-	-	-	-	-	-	-
10	⅜"		8	7,75	25	63,5	-	-	-	-	-	-	10,00	34	88,9	-	-	-	-	-	-	-	-	-	-	-	-	-	-	-
15	½"		8	9,40	25	63,5	-	-	-	9,40	25	108	-	-	-	-	-	-	-	-	-	-	-	-	-	-	-	-	-	-
10	⅜"	10	12,5	-	-	-	14,00	25,4	108	-	-	-	10,00	34	108	-	-	-	-	-	-	-	-	-	-	-	-	-	-	-
15	½"		12,5	9,40	25	88,9	18,10	50,5	108	9,40	25	108	16,00	34	108	-	-	-	-	-	-	-	-	-	-	-	-	-	-	-
20	¾"		12,5	15,75	25	101,6	-	-	-	15,75	25	117	-	-	-	-	-	-	-	-	-	-	-	-	-	-	-	-	-	-
15	½"	25	19	9,40	25	101,6	18,10	50,5	108	9,40	25	108	16,00	34	108	-	-	-	-	-	-	-	-	-	-	-	-	-	-	-
20	¾"		19	15,75	25	101,6	23,70	50,5	117	15,75	25	117	20,00	34	117	-	-	-	-	-	-	-	-	-	-	-	-	-	-	-
25	1"		19	22,10	50,5	114,3	29,70	50,5	127	22,10	50,5	127	26,00	50,5	127	22,60	50,5	127	23,00	50,5	127	23,00	50,5	127	-	-	-	-	-	-
32	1 ¼"	40	26	-	-	-	38,40	64	146	-	-	-	32,00	50,5	146	31,30	50,5	146	29,40	50,5	146	29,40	50,5	146	-	-	-	-	-	-
40	1 ½"		26	34,80	50,5	139,7	44,30	64	159	34,80	50,5	159	38,00	50,5	159	35,60	50,5	159	35,70	50,5	159	35,70	50,5	159	-	-	-	-	-	-
50	2"		32	47,50	64	158,75	56,30	77,5	190	47,50	64	190	50,00	64	190	48,60	64	190	47,80	64	190	47,80	64	190	-	-	-	-	-	-
65	2 ½"	80	62	60,20	77,5	193,68	72,10	91	216	60,20	77,5	216	66,00	91	216	60,30	77,5	216	59,50	77,5	216	59,50	77,5	216	-	-	-	-	-	-
80	3"		62	72,90	91	222,25	84,30	106	254	72,90	91	254	81,00	106	254	72,90	91	254	72,30	91	254	72,30	91	254	-	-	-	-	-	-
100	4"		76	97,38	119	292,1	109,70	144,5	305	97,38	119	305	100,00	119	305	97,60	119	305	97,60	119	305	97,60	119	305	-	-	-	-	-	-

Dimensions in mm

MG = diaphragm size

* Code 88 and Code 80 have different lengths

Materials and certificates

Tank valves are produced at GEMÜ from block material. The most frequently used material is 1.4435 (316L) and 1.4435 BN2 with a ferrite content of < 0.5%. In addition, other common alloys such as 1.4539, Titanium, Alloy 59, Hastelloy C and special alloys are available by agreement

The table below provides an overview of the possible certificates which are generally available. The type of certificate and its content must be specified exactly before ordering to be able to provide the required documents. Later requests of certificates may not be possible or possible only under certain conditions.

Our specialists are happy to answer any questions you might have.

Type	Designation of the test certificate in accordance with EN 10204	Content of the certificate	Confirmation of the certificate by
2.1	Certificate of compliance with the order	Confirmation of compliance with the order	the manufacturer
2.2	Test report	Confirmation of compliance with the order with specification of results of non-specific testing	the manufacturer
3.1	Inspection certificate 3.1	Confirmation of compliance with the order with specification of results of specific testing	the manufacturer acceptance officer independent of the production division
3.2	Inspection certificate 3.2	Confirmation of compliance with the order with specification of results of specific testing	the manufacturer acceptance officer independent of the production division and the acceptance officer commissioned by the purchaser or the acceptance officer named in the official regulations



Certificates and approvals

There is no universal diaphragm for all applications, hence we use different rubber mixtures and materials for our diaphragms. The different diaphragms have been certified according to their main fields of application so that we can certify conformity with the international regulations and rules below.

FDA (U.S. Food and Drug Administration)

FDA Extraction according to 21CFR 177.2600 for elastomers and 21CFR 177.1550 for PTFE.

USP (United States Pharmacopeia)

An independent institution has tested our diaphragms according to the regulations of the USP Class VI Chapter 87 In-Vitro and Chapter 88 In-Vivo. Furthermore, our diaphragms are free from animal ingredients.

RoHS

GEMÜ diaphragms comply with the RoHS Directive 2011/65 EC and the WEEE Directive 2002/96 EC.

Pressure Equipment Directive


As all diaphragm valves are pressure bearing components and as the diaphragm is the central sealing element in addition to the valve body, all diaphragms also comply with the European Pressure Equipment Directive 2014/68/EU Art. 3 § 3. If no original GEMÜ shut-off diaphragms are installed, GEMÜ cannot accept any responsibility.

EAC

Certification of conformity with the standards, quality and safety standards of the customs union between Russia, Kazakhstan, Belarus, Armenia and Kyrgyzstan.

EHEDG

Hygiene standards in the foodstuff and beverage industry are continuously increasing and are approaching those of the pharmaceutical industry in sensitive areas. For this reason, the EHEDG was established some years ago (European Hygienic Engineering and Design Group). GEMÜ was the first diaphragm valve manufacturer worldwide to be granted certification for its diaphragm valve system. The examination took place in 2002.

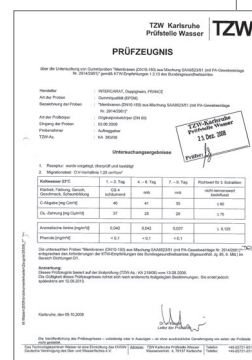

The original GEMÜ seal system

As a recognised diaphragm valve specialist, GEMÜ are familiar with almost all industrial sectors and applications. We are the leading supplier of stainless steel valves for aseptic and sterile applications in the pharmaceutical industry, biotechnology industry, as well as the foodstuff and beverage industries. As well as this, our valves also stand for reliability and a high standard of quality in the chemical and processing industries. The diaphragm, a central sealing element in the piping system is of major importance. Only the diaphragm and the valve body are in contact with the medium. At the same time, they also guarantee external hermetic sealing of the pipeline.


The system is more than the sum of the individual parts

The outstanding characteristics of the diaphragm valve are the result of the perfect interaction of tuned components. These are the valve body, the shut-off diaphragm, the diaphragm fixing, the compressor as well as the actuator. Our many years of experience and intensive dialogue with plant operators has enabled us to continue optimising the diaphragm valve design and its individual components.

Diaphragm and valve body are „inseparable“

GEMÜ valve bodies have a raised circular sealing bead on the inside diameter, in contrast to the valve bodies of other manufacturers. This results in a defined sealing edge. This measure reduces the ring-shaped gap between diaphragm and valve body in the external sealing area. This special feature makes GEMÜ diaphragm valves suitable for sterile applications. We also consider this crucial design and functional characteristic, which was developed by GEMÜ more than three decades ago and constantly refined, during the development of our diaphragms. Only this ensures that our customers can rely on the valve as a complete unit.

GEMÜ diaphragms have been developed, tested, and approved for applications with GEMÜ valve bodies. Therefore GEMÜ do not recommend or guarantee the use of other manufacturers diaphragms with GEMÜ valve bodies. We shall not accept any liability resulting from the use of diaphragms of other manufacturers inside GEMÜ diaphragm valves.


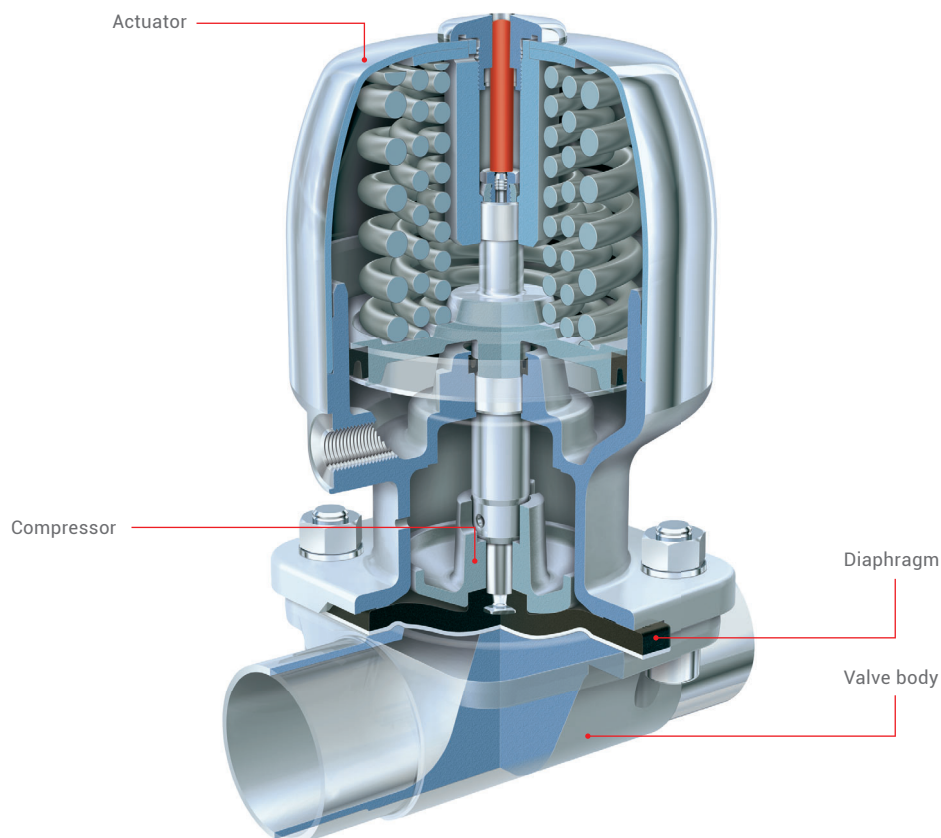


Flexible diaphragm fixing




The diaphragm is uniformly fixed in the compressor by means of a threaded pin. The only exception is the smallest diaphragm size (diaphragm size 8), which is pushed in with a rubber pin. The uniform fixing method applies both to soft elastomer and PTFE diaphragms. The largest advantage of fixing by means of a threaded pin, e.g. in comparison to a bayonet fitting, is the even transfer of forces onto the large area of the flanks of the screw thread. This prevents damage to the mechanical connection between compressor and diaphragm especially under vacuum operating conditions. The uniform fixing of elastomer and PTFE diaphragms enables subsequent replacement of the diaphragm while using the same actuator.




Diaphragm size 8

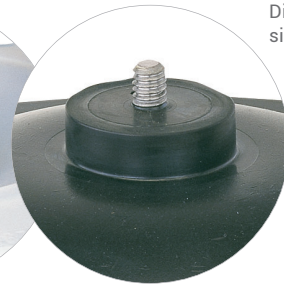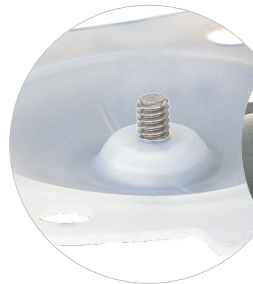

Diaphragm sizes 10 - 100

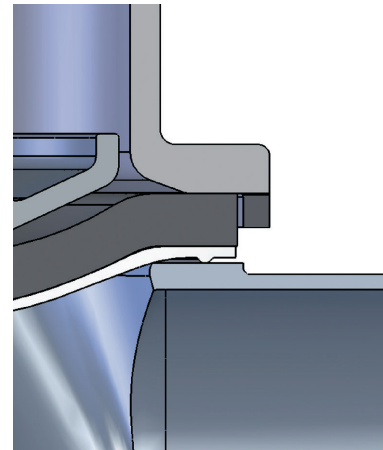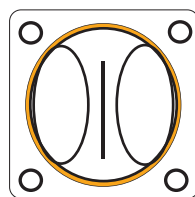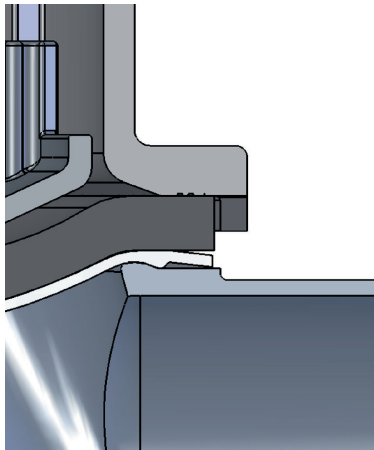

GEMÜ seal system

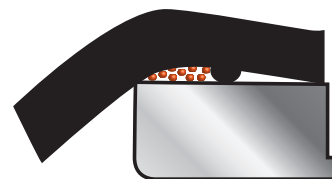

Conventional seal systems


GEMÜ diaphragms

Selection

			Temperature range [°C]			
Diaphragm	Material/Design	Diaphragm size	Liquid media			
			Min.	Max.	Sterilisation ¹	Code
EPDM	Ethylene-propylene-diene rubber	8 - 100	-10	100	max. 150 °C ² max. 60 min. per cycle	13/3A
EPDM	Ethylene-propylene-diene rubber	8 - 100	-10	100	max. 150 °C ² max. 180 min. per cycle	17
PTFE/EPDM	Fully laminated PTFE diaphragm with EPDM back	8, 10, 100	-10	100	max. 150 °C ² , no time limit per cycle	52/5A
PTFE/EPDM	Convex two-piece PTFE diaphragm with loose EPDM back	25, 40, 50, 80	-10	100	max. 150 °C ² , no time limit per cycle	5E

¹ The sterilisation temperature is valid for steam (saturated steam) or superheated water.

² If the sterilisation temperatures listed above are applied to the EPDM diaphragms for longer periods of time, the service life of the diaphragms will be reduced. In these cases, maintenance cycles must be adapted accordingly. This also applies to PTFE diaphragms exposed to high temperature fluctuations. PTFE diaphragms can also be used as moisture barriers; however, this will reduce their service life. The maintenance cycles must be adapted accordingly. GEMÜ 555 and 505 globe valves are particularly suitable for use in the area of steam generation and distribution. The following valve arrangement for interfaces between steam pipes and process pipes has proven itself over time: A globe valve for shutting off steam pipes and a diaphragm valve as an interface to the process pipes.

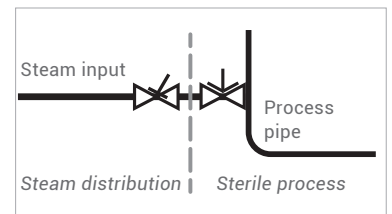

Each application must be analysed before the selection of the diaphragm material. Since the most varied operating conditions often prevail within a plant at different locations, it can be necessary to use different valves and materials. In particular, the chemical characteristics and the temperature of the working media often lead to different interactions. The suitability of the materials used must therefore always be examined individually with regard to the current resistance list or checked by an authorised specialist. Only this procedure guarantees that the application will operate safely and economically for a longer period. Diaphragms are wearing parts. They need to be regularly inspected and replaced otherwise malfunctions can occur, possibly resulting in hazardous situations. Please note: The maintenance intervals for inspecting and replacing diaphragms are application-dependent. In order to determine a suitable maintenance interval, the maintenance history and the stresses placed on the parts due to frequent sterilisation or frequent cycle duties must be taken into account.

Note

Since plastics and elastomers are subject to natural aging, we recommend observing the GEMÜ storage conditions for shut-off diaphragms. You thereby guarantee maximum storage and service life of the diaphragms. The temperatures specified above are merely the permissible temperature ranges for the respective diaphragm. The permissible temperature ranges of the valve must always be taken into account for the overall valve design. These can be found in the respective datasheets. The temperature values are indicated irrespective of operating pressure and diaphragm size and apply to water and/or inert gases. When using water vapour or saturated steam, observe the steam pressure diagram.

Certificates and approvals





FDA compliant	USP Class VI	EHEDG	TA Luft (German Clean Air Act)	O ₂ BAM	Compatibility with media	Special features
•	•	•	•	•	Very good all-round elastomer, resistant to many acidic and alkaline media, demineralised and deionised hot water, inert and many other industrial gases.	Suitable for vacuum, low gas permeability, applicable for steam sterilisation.
•	•	•	•	•		Compound and construction of the diaphragm have been specially optimised for steam applications, clearly improved service life.
•	•	•	•	•	Resistant to nearly all chemicals, such as strong acids, alkalis and salts, also at high temperatures, steam, WFI as well as pharmaceuticals. Good resistance to solvents, chlorine, and aromatic hydrocarbons.	Fully laminated diaphragm, can be used in steam. Low gas permeability.
•	•	•	•	•		Convex two-piece diaphragm with loose PTFE face for higher switching cycles, can be used for permanent steam application. Special compounding and production by GEMÜ. Special seal contour for external sealing on the bottom of the diaphragm. Low gas permeability.

EPDM diaphragm dimensions [mm]

MG*	DN	NPS	A	B	C	D	ød	e	h	W	α	β	γ	Y	Number of bolt holes
8	4 - 15	¼" - ½"	22	22	31,5	31,5	4,5	4	5,6	-	-	-	-	-	4
10	10 - 20	⅜" - ¾"	39	44	48	53	5,2	5	9	M4	-	-	-	-	4
25	15 - 25	½" - 1"	54	46	71,7	66,7	9	6	8	¼"	-	-	-	-	4
40	32 - 40	1¼" - 1½"	70	65	100	90	11,5	7	8	¼"	-	-	-	-	4
50	50	2"	82	78	124	106	13	7	7	¼"	-	-	-	-	4
80	80	3"	127	114	186	156	18	9	8	5/16"	-	-	-	-	4
100	100	4"	194	-	228	-	13	10	9	5/16"	28°	42°	40°	-	8

* Diaphragm size

The thread of the diaphragm pin „W" corresponds to Whitworth standard.

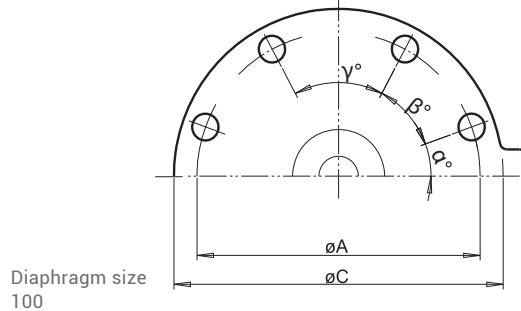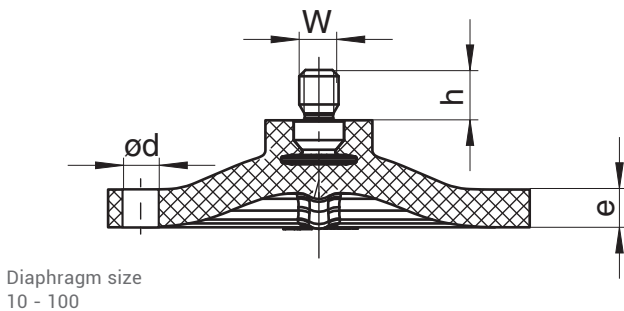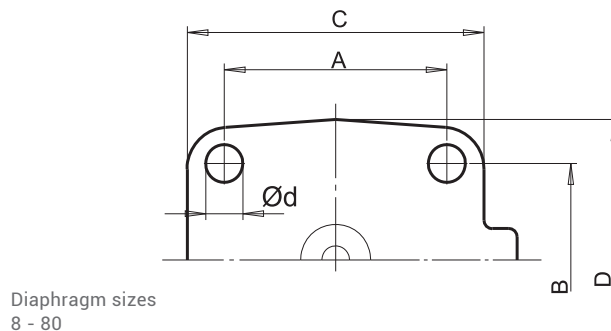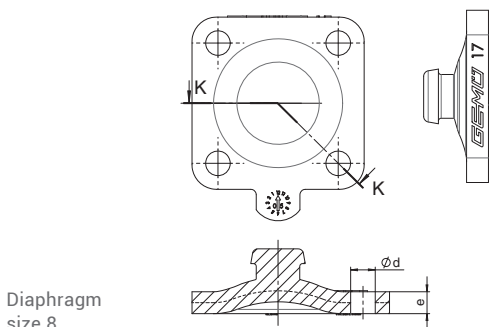

GEMÜ B600 tank valves

Selection of operators

Manually operated



Type	GEMÜ 9601	GEMÜ 9602	GEMÜ 9612	GEMÜ 9673	GEMÜ 9653	GEMÜ 9654
Material	Stainless steel, plastic handwheel, with optical position indicator and seal adjuster	Stainless steel, with optical position indicator and seal adjuster	Stainless steel, plastic handwheel, with optical position indicator and seal adjuster	Stainless steel, plastic handwheel, with optical position indicator and seal adjuster	Stainless steel, plastic handwheel, with optical position indicator, stroke limiter/seal adjuster, lockable, optional: electrical position indicator	Stainless steel, with optical position indicator, stroke limiter/seal adjuster, lockable, optional: electrical position indicator
Autoclavable	●	●	●	●	●	●
Operating temperature*	-10 to 150 °C	-10 to 150 °C	-10 to 150 °C	-10 to 150 °C	-10 to 150 °C	-10 to 150 °C
Operating pressure*	0 to 10 bar	0 to 10 bar	0 to 10 bar	0 to 10 bar	0 to 10 bar	0 to 10 bar
DN	4 to 15	4 to 15	10 to 20	15 to 50	10 to 100	4 to 100
Diaphragm size 8	●	●	-	-	-	●
Diaphragm size 10	-	-	●	-	●	●
Diaphragm size 25	-	-	-	●	●	●
Diaphragm size 40	-	-	-	●	●	●
Diaphragm size 50	-	-	-	●	●	●
Diaphragm size 80	-	-	-	-	●	●
Diaphragm size 100	-	-	-	-	●	●

* dependent on diaphragm material, see technical datasheet

Elastomer diaphragms

EPDM



PTFE diaphragms

PTFE/EPDM, PTFE/FPM




Pneumatically operated

							
GEMÜ 9605	GEMÜ 9625	GEMÜ 9687	GEMÜ 9650	GEMÜ 9650TL	GEMÜ 9651	GEMÜ 9658/9688	GEMÜ 9660
Plastic, with stainless steel distance piece, optical position indicator	Plastic, with stainless steel distance piece, optical position indicator	Plastic, with stainless steel distance piece	Stainless steel, with optical position indicator, optionally autoclavable	Safety valve, stainless steel, mounting facility for proximity switches	Stainless steel, with integrated automation module	Two stage actuator, stainless steel	Filling valve, stainless steel with optical position indicator
-	-	-	● (DN 4 to 25)	-	-	-	-
-10 to 150 °C	-10 to 150 °C	-10 to 150 °C	-10 to 150 °C	-10 to 150 °C	-10 to 150 °C	-10 to 150 °C	-10 to 150 °C
0 to 8 bar	0 to 6 bar	0 to 10 bar	0 to 10 bar	0 to 8 bar	0 to 10 bar	0 to 10 bar	0 to 5 bar
4 to 15	10 to 20	10 to 100	4 to 100	4 to 25	4 to 25	10 to 50	4 to 25
●	-	-	●	●	●	●	●
-	●	●	●	●	●	●	●
-	-	●	●	●	●	●	●
-	-	●	●	-	-	●	-
-	-	●	●	-	-	●	-
-	-	●	●	-	-	-	-
-	-	●	●	-	-	-	-

Other valve body versions



2/2-way body investment casting
2/2-way version to all international standard butt weld connections

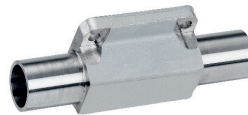

2/2-way body forged version
2/2-way version to all international standard butt weld connections

Other connections



Clamps
to all common standards



Aseptic clamps
to all common standards



Aseptic unions
to all common standards



Aseptic flanges
to all common standards

Other versions, accessories as well as motorized operators available.

GEMÜ B600 tank valves

Selection of operators

Motorized



Typ	GEMÜ 9618	GEMÜ 9698
Material	Plastic, with stainless steel distance piece, optical position indicator	Plastic, with stainless steel distance piece, optical position indicator and manual override
Autoclavable	-	-
Operating temperature*	0 to 130 °C (ohne Zwischenstück 15 to 50 °C)	-10 to 150 °C
Operating pressure*	0 to 6 bar	0 to 6 bar
DN	4 - 15	15 - 50
Anschlussspannung	24 VAC, 120 VAC, 230 VAC, 50/60Hz	24 VAC, 120 VAC, 230 VAC, 50/60Hz
Diaphragm size 8	●	-
Diaphragm size 10	●	-
Diaphragm size 25	-	●
Diaphragm size 40	-	●
Diaphragm size 50	-	●
Diaphragm size 80	-	-
Diaphragm size 100	-	-

* dependent on diaphragm material, see technical datasheet

Positioners and process controllers

Overview

GEMÜ pneumatic actuators can usually be paired with our electro-pneumatic positioners and process controllers.

In addition to the process parameters and the control system for which a positioner must be suitable, other technical functions and properties also play an important part in the selection of the right positioner.

To make your choice easier, we have placed the four GEMÜ positioners in a comparison based on important features.



Function / Features		GEMÜ 1434 µPos	GEMÜ 1435 ePos	GEMÜ 1436 cPos	GEMÜ 1436 cPos eco
Controller type	Positioner	•	•	•	•
	Process controller			•	
Operation	Local display / keypad		•	•	
	Status display	•	•	•	•
	Web browser user			•	
	Fieldbus option (Profibus DP, Device Net)			•	
Housing	Plastic	•		•	•
	Aluminium		•		
Functions	Automatic initialisation (speed ^{-AP})	•	•	•	•
	Alarm / error outputs		•	•	
	Min/max positions adjustable		•	•	
Mounting	Direct mounting to linear actuators	•	•	•	•
	Remote mounting to linear actuators	•	•	•	•
	Direct mounting to quarter turn actuators		•	•	•
	Remote mounting to quarter turn actuators		•	•	•
Control function of valve actuator	Control function 1, normally closed (NC)	•	•	•	•
	Control function 2, normally open (NO)	•	•	•	•
	Control function 3, double acting (DA)		•	•	
Air output		15 NI/min.	50 NI/min. 90 NI/min.	150 NI/min. 200 NI/min. 300 NI/min.	150 NI/min. 200 NI/min.

Combi switchboxes and electrical position indicators for pneumatically operated linear valves

Our devices detect the valve stroke in any installation position without play and are tension-free. The sensor base of the GEMÜ 1234, 1235 and 4242 series is positively connected to the valve spindle by means of a preloaded spring so that possible tangential forces of the valve actuator do not negatively affect the position indicator. The position indicators can be quickly and easily assembled and are safe and uncomplicated to handle.

They can be adapted to pneumatic actuators of GEMÜ globe and diaphragm valves. Due to its special design, the GEMÜ 9415 quarter turn actuator may likewise be combined with this type of position indicator.



Device type	Combi switchbox Electrical position indicators										
	4242	1215	1230	1231	1232	1201	1211	1214	1234	1235/ 1236	1242
Valve stroke (in mm)	2 - 75		2 - 20	2 - 20	2 - 20	2 - 70	2 - 70	2 - 70	1 - 10	2 - 30 4 - 50 5 - 75	2 - 46
Electrical connection	M12	¹⁾	¹⁾	²⁾	¹⁾	²⁾	²⁾	¹⁾	M12	M12	M12
Programmable	•								•	•	•
With integrated pilot valve	•										
Ex design		• ³⁾		•			•				
NEC 500 version	•										•
Fieldbus interface	•								•		•
Mechanically adjustable switches			•	•	•	•	•	•			
Mechanically adjustable (microswitches)			•			•					
UL version	•		•								•
SIL version										•	
Optical position indicator (LEDs)	•		•		•			•	•	•	•
Optical position indicator (mechanical)	• ⁴⁾	•									
Feedback (OPEN and CLOSED)	•		•	•	•	•	•	•	•	•	•
Feedback (OPEN)		•									

¹⁾ cable gland, M12 (optional)

²⁾ cable gland

³⁾ not with optional M12 plug

⁴⁾ for size 2 only

GEMÜ B600 tank valves

Specification

The enormous consequential costs which may be incurred as a result of errors in the planning of production plant more than justify increased planning efforts. Delays and extra costs in validation, late commissioning, contaminated batches, later modifications to the plant are just a few of the points in favour of precise planning.

Good planning of the valve designs begins in the project stage. The implementation of complex process sequences demands a wide variety of compact valve designs. GEMÜ places great emphasis on ensuring you get the optimum tank valve for your specific application and not any standard solution from the catalogue. It is therefore important that you give us all the relevant information in the project phase. The GEMÜ specification sheet is a great help here.

Please use only one form for each tank valve and proceed as follows to fill in the specification:

1. Enter the operating conditions and desired materials.
2. Please state what functions the tank valve should fulfil. Draw a pictogram or functional diagram and insert it in the specification. You can of course use the examples shown in this brochure as a guide.
3. Label all connection spigots starting with S1, all valve seats starting with V1.
4. Assign the necessary features to every connection in the table and add explanatory remarks where necessary.
5. Specify the necessary operator and type as well as control function for every connection.
6. For extra remarks and descriptions, you can use an additional sheet.



Specification of B600 tank valves

Please complete this form and return it to your nearest GEMÜ office or to the address listed below!

Operating pressure: bar

Medium temperature: °C

Material of tank valve body:

1.4435 ☐

1.4435 BN 2 ($\Delta Fe < 0,5\%$) ☐

1.4539 ☐

Other ☐

Tests/Certificates:

AD 2000 W2 (Standard) ☒

Other ☐

Diaphragm material:

EPDM Code

PTFE Code

Other

Surface of tank valve body:

1502 (Ra) $\leq 0.8 \mu m$ ☐

1503 (Ra) $\leq 0.8 \mu m$ electropolished ☐

1507 (Ra) $\leq 0.6 \mu m$ ☐

1508 (Ra) $\leq 0.6 \mu m$ electropolished ☐

1536 (Ra) $\leq 0.4 \mu m$ ☐

1537 (Ra) $\leq 0.4 \mu m$ electropolished ☐


1527 (Ra) $\leq 0.25 \mu m$ ☐

1516 (Ra) $\leq 0.25 \mu m$ electropolished ☐

Other ☐

Quantity:

Example: B600 03-02.A

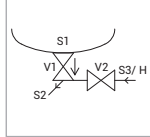


Please draw functional diagram.

Important: Please observe correspondence of table and functional diagram.

Please specify design (e.g. B600 03-02.A):

Tank radius R= mm

Welding neck thickness SP = mm
(Standard 6 mm)

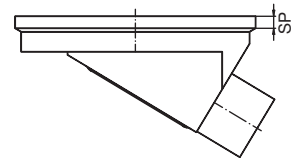

Draining direction:

Spigot: S1, S2, ...

Valve seat:

Preferred installation position: Horizontal/Vertical

Intersection:

Flow direction (medium):

Spigot	Pipe connection				Operator			Other
Spigot no.	DN	Code	$\phi d(a)[mm]$	s[mm]	Operator type	Control function	Operator size	Comment/accessories
S1	Welding diameter dependent on type and diaphragm size							
S2					V1			
S3					V2			
S4					V3			
S5					V4			

The technical details of each enquiry must be checked by GEMÜ.

Contact (GEMÜ):

Customer:

Department:

Address:

Phone: E-mail:

Please do not write here!

K-No.:

P600:

B600:

X:

Worldwide presence

AUSTRALIA

GEMÜ Australia Pty. Ltd
Unit 4 - 8/10 Yandina Road
West Gosford, NSW 2250
Phone: +61-2-43 23 44 93
Fax: +61-2-43 23 44 96
mail@gemu.com.au

AUSTRIA

GEMÜ GmbH
Europaring F15 401
2345 Brunn am Gebirge
Phone: +43 2236 30 43 45-0
Fax: +43 2236 30 43 45-31
info@gemue.at

BELGIUM

GEMÜ Valves bvba/sprl
Koning Albert 1 laan, 64
1780 Wemmel
Phone: +32 2 702 09 00
Fax: +32 2 705 55 03
info@gemue.be

BRAZIL / SOUTH AMERICA

GEMÜ Indústria de Produtos
Plásticos e Metalúrgicos Ltda.
Rue Marechal Hermes, 1141
83.065-000 São José dos Pinhais
Paraná
Phone: +55-41-33 82 24 25
Fax: +55-41-33 82 35 31
gemu@gemue.com.br

CANADA

GEMÜ Valves Canada Inc.
2572 Daniel-Johnson Boulevard
Laval, Quebec
H7T 2R8
Phone: +1-450-902-2690
Fax: +1-404-3 44 4003
info@gemu.com

CHINA

GEMÜ Valves (China) Co., Ltd
No.518, North Hengshahe Road
Minhang District, 201108
Shanghai
Phone: +86-21-2409 9878
info@gemue.com.cn

DENMARK

GEMÜ ApS
Industriparken 16-18
2750 Ballerup
Phone: +45 70 222 516
info@gemue.dk

FRANCE

GEMÜ S.A.S
1 Rue Jean Bugatti
CS 99308 Duppigheim
67129 Molsheim Cedex
Phone: +33-3 88 48 21 00
info@gemu.fr

INTERCARAT
1 Rue Jean Bugatti
CS 99308 Duppigheim
67129 Molsheim Cedex
Phone: +33-3 88 48 21 20
sales@intercarat.com

GERMANY

GEMÜ Gebr. Müller GmbH & Co. KG
Fritz-Müller-Straße 6 - 8
74653 Ingelfingen-Criesbach
Postfach 30
74665 Ingelfingen-Criesbach
Phone: +49 (0)7940-12 30
Fax: +49 (0)7940-12 31 92
(Domestic)
Fax: +49 (0)7940-12 32 24 (Export)
info@gemue.de

Inevvo solutions GmbH & Co. KG
Fritz-Müller-Platz 1
74676 Niedernhall-Waldzimmern
Phone: +49 (0)7940-12 38 681
info@inevvo-solutions.com

GREAT BRITAIN

GEMÜ Valves Ltd.
10 Olympic Way
Birchwood, Warrington
WA2 0YL
Phone: +44-19 25-82 40 44
Fax: +44-19 25-82 80 02
info@gemu.co.uk

HONG KONG

GEMÜ (Hong Kong) Co., Ltd.
Room 2015, Tower B,
Regent Centre,
70 TA Chuen Ping Street
Kwai Chung, N.T., Hong Kong
P.R. China
Phone: +852 6873 8280
Fax: +852 6873 8280
info@gemue.com.cn

INDIA

GEMÜ Branch Office India
Room Number 135,
1st Floor, 101-104, B-Wing,
GCP Business Centre
Opp. Memnagar Fire Station
Vijay Cross Road
Ahmedabad - 380 014
Phone: +91-79-6134 4423
Fax: +91-79-25450439
sales@gemu.in

INDONESIA

GEMÜ Valves Pte Ltd
(Indonesia Representative Office)
Rukan Mangga Dua Square
Block F17, 2nd Floor
Jl. Gunung Sahari Raya No. 1
Jakarta Utara 14420
Indonesia
Phone: +62 (21) - 6231 0035
Fax: +62 (21) - 2907 4643
info@gemu.co.id

IRELAND

GEMÜ Ireland Ltd
15 Eastgate Drive
Eastgate Business Park
Little Island
Co. Cork
Phone: +353 (0)21 4232023
Fax: +353 (0)21 4232024
info@gemu.ie

ITALY

GEMÜ S.r.l.
Via Giovanni Gentile, 3
20157 Milano
Phone: +39-02-40044080
Fax: +39-02-40044081
info@gemue.it

JAPAN

GEMÜ Japan Co., Ltd.
2-5-6, Aoi, Higashi-ku,
Nagoya, Aichi, 461-0004
Phone: +81-52-936-2311
Fax: +81-52-936-2312
info@gemu.jp

MALAYSIA

GEMÜ Valves PTE. LTD.
D-8-09, Block D, No. 2A
Jalan PUJ 1A/7A
Oasis Square, Oasis Damansara
47301 Petaling Jaya
Selangor Darul Ehsan
Phone: +(603)- 7832 7640
Fax: +(603)- 7832 7649
info@gemu.com.sg

MEXICO

GEMÜ Valvulas S.A. de C.V.
German Centre,
Av. Santa Fe No. 170 - OF. 5-1-05
Col. Lomas de Santa Fe,
Del. Alvaro Obregon
01210 Mexico, D.F.
Phone: +52 55 7090 4161
+52 55 7090 4179

RUSSIA

OOO „GEMÜ GmbH“
Uliza Shipilovskaya, 28A
115563, Moskau
Phone: +7(495) 662-58-35
Fax: +7(495) 662-58-35
info@gemue.ru

SINGAPORE

GEMÜ Valves PTE. LTD.
25 International Business Park
German Centre #03-73/75
Singapore 609916
Phone: +65-65 62 76 40
Fax: +65-65 62 76 49
info@gemu.com.sg

SOUTH AFRICA

GEMÜ Valves Africa Pty. Ltd
Cnr Olympic Duel Avenue
And Angus Crescent,
Northlands Business Park
(Stand 379),
New Market Road
Randburg
Phone: +27 11 462 7795
Fax: +27 11 462 4226
info@gemue.co.za

SWEDEN

GEMÜ Armatur AB
Heljesvägen 8
437 36 Lindome
Phone: +46-31-996500
order@gemu.se

SWITZERLAND

GEMÜ GmbH
Seetalstr. 210
6032 Emmen
Phone: +41-41-7 99 05 05
Fax: +41-41-7 99 05 85
info@gemue.ch

GEMÜ Vertriebs AG
Lettenstrasse 3
6343 Rotkreuz
Phone: +41-41-7 99 05 55
Fax: +41-41-7 99 05 85
vertriebsag@gemue.ch

TAIWAN

GEMÜ Taiwan Ltd.
9F.-5, No.8, Ziqiang S. Rd.
Zhubei City
Hsinchu County 302,
Taiwan (R.O.C.)
Phone: +886-3-550-7265
Fax: +886-3-550-7201
office@gemue.tw

UNITED STATES

GEMÜ Valves Inc.
3800 Camp Creek Parkway
Suite 120, Building 2600
Atlanta, Georgia 30331
Phone: +1-678-5 53 34 00
Fax: +1-404-3 44 93 50
info@gemu.com

In addition to these
subsidiaries, GEMÜ has a
global partner network.

Contact details:
[www.gemu-group.com/
en_GB/kontakte/](http://www.gemu-group.com/en_GB/kontakte/)



 GEMÜ manufacturing site

 GEMÜ subsidiary

